

# memmert

UN<sub>xx</sub>plus | UF<sub>xx</sub>plus  
IN<sub>xx</sub>plus | IF<sub>xx</sub>plus  
UN<sub>xx</sub>mplus | UF<sub>xx</sub>mplus  
IN<sub>xx</sub>mplus | IF<sub>xx</sub>mplus



## OPERATING INSTRUCTIONS

UNIVERSAL OVEN | U  
INCUBATOR | I

MADE IN GERMANY.

[www.memmert.com](http://www.memmert.com)

## Manufacturer and customer service

Memmert GmbH + Co. KG  
Willi Memmert Straße 90-96  
D-91186 Büchenbach  
Deutschland

Phone: +49 (0)9122 925-0  
Fax: +49 (0)9122 14585  
E-mail: [sales@memmert.com](mailto:sales@memmert.com)  
Internet: [www.memmert.com](http://www.memmert.com)

### Customer service:

Service hotline: +49 (0)9171 9792 911  
Service fax: +49 (0)9171 9792 979  
E-mail: [service@memmert.com](mailto:service@memmert.com)

When contacting customer service, always quote the product serial number on the nameplate (see page 13).

### Shipping address for repairs:

Memmert GmbH + Co. KG  
Kundenservice  
Willi-Memmert-Str. 90-96  
DE-91186 Büchenbach  
Germany

Please contact our customer service before sending appliances for repair or before returning equipment, otherwise, we have to refuse acceptance of the shipment.

© 2020 MEMMERT GmbH + Co. KG

D33355 | Date 01/2020

We reserve the right to make changes

## About this manual

---

### Purpose and target group

This manual describes the assembly, function, transport and operation of the universal ovens UN<sub>x</sub>plus/UF<sub>x</sub>plus/UN<sub>x</sub>mplus/UF<sub>x</sub>mplus and the incubators IN<sub>x</sub>plus/IF<sub>x</sub>plus/IN<sub>x</sub>mplus/IF<sub>x</sub>mplus. It is intended for use by trained personnel of the owner, who have the task of operating and/or maintaining the respective appliance.

If you are asked to work on the appliance, read this manual carefully before starting. Familiarise yourself with the safety regulations. Only perform work that is described in this manual. If there is something you do not understand, or certain information is missing, ask your superior or contact the manufacturer. Do not do anything without authorisation.

### Versions

The appliances are available in different configurations and sizes. If specific equipment features or functions are available only for certain configurations, this is indicated at the relevant points in this manual.

The functions described in this manual refer to the latest firmware version.

Due to individual configurations and sizes, illustrations in this manual may be slightly different from the actual appearance. Function and operation are identical.

### Other documents that have to be observed:

- ▶ For operation of the appliance with MEMMERT AtmoCONTROL, observe the respective software manual. To open the AtmoCONTROL software manual, click on "Help" in the AtmoCONTROL menu bar.
- ▶ For service and repair (see page 56), please refer to the separate service manual

### Storage and forwarding

This instruction manual belongs with the appliance and should always be stored where persons working on the appliance have access to it. It is the responsibility of the owner to ensure that persons who are working or will work on the appliance are informed as to the whereabouts of this instruction manual. We recommend that it is always stored in a protected location close to the appliance. Make sure that the instruction manual is not damaged by heat or humidity. If the appliance is sold on or transported and then set up again at a different location, the operating instructions must go with it.

You will find the current version of our operating manual as pdf file if you go to [www.memmert.com/de/service/downloads/bedienungsanleitung/](http://www.memmert.com/de/service/downloads/bedienungsanleitung/).

# Contents

---

<b>1.</b>	<b>For your Safety</b>	<b>6</b>
1.1	Terms and signs used.....	6
1.2	Product safety and dangers .....	7
1.3	Requirements of the operating personnel .....	7
1.4	Responsibility of the owner.....	8
1.5	Changes and alterations.....	8
1.6	Behaviour in case of malfunctions and irregularities.....	8
1.7	Switching off the appliance in an emergency .....	8
<b>2.</b>	<b>Construction and description</b>	<b>9</b>
2.1	Construction .....	9
2.2	Function.....	11
2.3	Material.....	11
2.4	Electrical equipment.....	11
2.5	Connections and interfaces.....	12
2.6	Designation (nameplate).....	12
2.7	Technical data .....	13
2.8	Declaration of conformity .....	15
2.9	Ambient conditions .....	15
2.10	Scope of delivery .....	16
2.11	Optional accessories.....	16
<b>3.</b>	<b>Delivery, transport and setting up</b>	<b>17</b>
3.1	For your Safety.....	17
3.2	Delivery .....	18
3.3	Transport.....	18
3.4	Unpacking .....	18
3.5	Storage after delivery .....	18
3.6	Setting up .....	19
<b>4.</b>	<b>Putting into operation</b>	<b>23</b>
4.1	Connecting the appliance .....	23
4.2	Switching on.....	24
<b>5.</b>	<b>Operation and control</b>	<b>25</b>
5.1	Operating personnel.....	25
5.2	Opening the door.....	25
5.3	Loading the appliance.....	26
5.4	Operating the appliance.....	26
5.5	Temperature monitoring .....	33
5.6	Graph.....	37
5.7	Ending operation.....	37



---

<b>6.</b>	<b>Malfunctions, warning and error messages</b>	<b>38</b>
6.1	Warning messages of the monitoring function.....	38
6.2	Malfunctions, operating problems and appliance errors .....	39
6.3	Power failure.....	41
<b>7.</b>	<b>Menu mode</b>	<b>42</b>
7.1	Overview.....	42
7.2	Basic operation in menu mode using the example of language selection.....	43
7.3	Setup.....	44
7.4	Date and Time .....	48
7.5	Calibration .....	50
7.6	Programme.....	52
7.7	Sound .....	53
7.8	Protocol .....	54
7.9	User ID .....	55
<b>8.</b>	<b>Maintenance and service</b>	<b>56</b>
8.1	Cleaning.....	56
8.2	Regular maintenance.....	56
8.3	Repairs and service .....	56
<b>9.</b>	<b>Storage and disposal</b>	<b>57</b>
9.1	Storage .....	57
9.2	Disposal .....	57

---

# 1. For your Safety

## 1.1 Terms and signs used

In this manual and on the appliance itself, certain common terms and signs are used to warn you of possible dangers or to give you hints that are important in avoiding injury or damage. Observe and follow these hints and regulations to avoid accidents and damage. These terms and signs are explained below.

### 1.1.1 Terms used

**"Warning"** is used whenever you or somebody else could be injured if you do not observe the accompanying safety regulation.

**"Caution"** is used for information that is important for avoiding damage.

### 1.1.2 Signs used

Warning signs (warning of a danger)					
Danger of electrocution	Danger of explosion	Dangerous gases / vapours	Danger of burns	Danger of toppling over	Hazard area! Observe the operating instructions
Prohibition signs (forbidding an action)					
Do not lift	Do not tilt	Do not enter			
Regulation signs (stipulating an action)					
Disconnect the mains plug	Wear gloves	Wear safety boots	Observe information in separate manual		
Other icons					
	Important or useful additional information				

## 1.2 Product safety and dangers

The appliances described in this manual are technically sophisticated, manufactured using high-quality materials and subject to many hours of testing in the factory. They contain the latest technology and comply with recognised technical safety regulations. However, there are still risks involved, even when the appliances are used as intended. These are described below.



### Warning!

After removing covers, live parts may be exposed. You may receive an electric shock if you touch these parts. Disconnect the mains plug before removing any covers. Only electrical technicians may work on the electrical equipment of the appliances.



### Warning!

When loading the appliance with an unsuitable load, poisonous or explosive vapours or gases may be produced. This could cause the appliance to explode, and people could be severely injured or poisoned. The appliance may only be loaded with materials/test objects which do not form any toxic or explosive vapours when heated up (see also page 10).



### Warning!

If the door is open while the appliance is in operation, the appliance may overheat and pose a fire hazard. Do not leave the door open during operation.



### Warning!

Depending on operation, the surfaces in the working chamber and the chamber load may still be very hot after the appliance is switched off. Touching these surfaces can cause burns. Wear heat-resistant protective gloves or wait until the appliance cools down.



### Warning!

In case of appliances of a certain size, you can get accidentally locked in, which is life-threatening. Do not climb into the appliance!

## 1.3 Requirements of the operating personnel

The appliance may only be operated and maintained by persons who are of legal age and have been instructed accordingly. Personnel who are to be trained, instructed or who are undergoing general training may only work with the appliance under the continuous supervision of an experienced person.

Repairs may only be performed by qualified electricians. The regulations in the separate service manual must be observed.

## 1.4 Responsibility of the owner

The owner of the appliance

- ▶ is responsible for the flawless condition of the appliance and for its proper operation in accordance with its intended use (see page 10);
- ▶ is responsible for ensuring that persons who are to operate or service the appliance are qualified to do this, have been instructed accordingly and are familiar with the operating instructions at hand;
- ▶ must know about the applicable guidelines, requirements and operational safety regulations, and train staff accordingly;
- ▶ is responsible for ensuring that unauthorised persons have no access to the appliance;
- ▶ is responsible for ensuring that the maintenance plan is adhered to and that maintenance work is carried out properly (see page 56);
- ▶ has to ensure that the appliance and its surroundings are kept clean and tidy, for example through corresponding instructions and inspections;
- ▶ is responsible for ensuring that personal protective clothing is worn by operating personnel, e.g. work clothes, safety shoes and protective gloves.

## 1.5 Changes and alterations

No unauthorised changes or alterations may be made to the appliance. No parts may be added or inserted which have not been approved by the manufacturer.

Unauthorised modifications or changes result in the CE declaration of conformity losing its validity and the appliance must no longer be operated.

The manufacturer is not liable for any damage, danger or injuries that result from unauthorised changes or alterations, or from non-observance of the regulations in this manual.

## 1.6 Behaviour in case of malfunctions and irregularities

The appliance may only be used in a flawless condition. If you as the operator notice irregularities, malfunctions or damage, immediately take the appliance out of service and inform your superior.


**i** You can find information on correcting malfunctions from page 38.

## 1.7 Switching off the appliance in an emergency

Push the On/Off switch on the control panel (Fig. 1) and disconnect power plug. This disconnects the appliance from the power supply at all poles.



**Warning!**  
Depending on operation, the surfaces in the working chamber and the chamber load may still be very hot after the appliance is switched off. Touching these surfaces can cause burns. Wear heat-resistant protective gloves or wait until the appliance cools down.

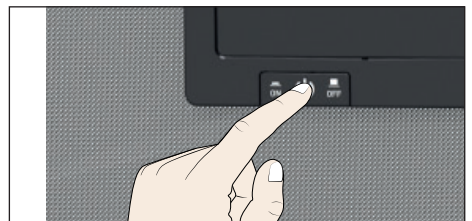

*Fig. 1*  
Switch off the appliance by pressing the On/Off switch

## 2. Construction and description

### 2.1 Construction



Fig. 2 Construction

- |  |                                      |
|--|--------------------------------------|
| 1 ControlCOCKPIT with capacitive function keys (see page 27) | 5 Working chamber                    |
| 2 On/Off switch (see page 24)                                | 6 Nameplate (covered, see page 13)   |
| 3 Working chamber fan (for UF.../IF... appliances only)      | 7 Door handle (see page 25)          |
| 4 Steel grid   | 8 Turn control with confirmation key |
|  | 9 USB interface (see page 12)        |

## 2.2 Intended use

Based on the standards and guidelines listed in the following, the products described in this manual have received a CE label from the company Memmert:



Product	Intended use	Directives
UN <sub>xx</sub> plus UF <sub>xx</sub> plus IN <sub>xx</sub> plus IF <sub>xx</sub> plus	<p>This appliance is exclusively intended for heating up non-explosive substances and objects. Any other use is improper and may result in hazards and damage.</p> <p>The appliance is not explosion-proof (does not comply with the German occupational health and safety regulation VBG 24). The appliance may only be loaded with materials and substances which cannot form any toxic or explosive vapours at the set temperature and which cannot explode, burst or ignite.</p> <p>The appliance may not be used for the drying, evaporation and burning-in of paints or similar materials, the solvents of which could form an explosive mixture when combined with air. If there is any doubt as to the composition of materials, they must not be loaded into the appliance. Potentially explosive gas-air mixtures must not form, neither in the working chamber nor in the direct vicinity of the appliance.</p>	<ul style="list-style-type: none"> <li>▶ Directive 2004/108/EC amended (Directive of the European Parliament and of the Council on the approximation of the laws of the Member States relating to electromagnetic compatibility). Standards complied with: DIN EN 61326:2004-05, EN 61326:1997, EN 61326/A1:1998, EN 61326/A2:2001, EN 61326/A2:2003</li> <li>▶ Directive 2006/95/EC amended (Directive of the European Parliament and of the Council on the harmonisation of the laws of Member States relating to electrical equipment designed for use within certain voltage limits). Standards complied with: DIN EN 61 010-1 (VDE 0411 part 1), DIN EN 61 010-2-010 (VDE 0411 part 2-010), EN 61 010-1:2001, EN 61 010-2-010</li> </ul>
Product	Intended use as a medical device	Directive
UN <sub>xx</sub> mplus UF <sub>xx</sub> mplus	The appliance may be used for heating fango, silicate and APS packs for physical therapy and keeping them warm.	Directive 93/42/EEC (Council Directive concerning medical devices)
IN <sub>xx</sub> mplus IF <sub>xx</sub> mplus	The appliance may be used for temperature control of rinsing and infusion solutions and contrast agents.	

## 2.3 Function

Appliances of the UN... and IN... type series feature natural circulation (convection). For the UF... and IF... type series, air is circulated with a fan at the working chamber rear panel (Fig. 3, No. 1). It increases the air flow and provides stronger horizontal forced air circulation than natural convection.

In both the convection and fan ventilated appliances, supply air (2) is preheated in a pre-heating chamber (3). Through the ventilation slits in the side panel of the working chamber, the pre-heated air is introduced into the interior of the chamber. The supply and exhaust air (5) volume (air change) is controlled by the air flap (4) on the rear panel of the appliance.

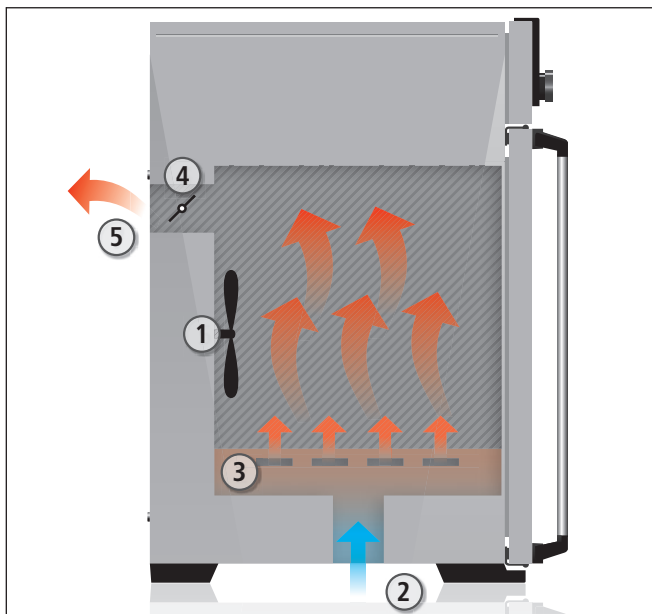


Fig. 3 Function

- 1 Fan
- 2 Fresh air
- 3 Pre-heating chamber
- 4 Air flap
- 5 Exhaust air

## 2.4 Material

For the outer housing, MEMMERT deploys stainless steel (Mat.No. 1.4016 – ASTM 430) and for the interior, stainless steel (Mat.No. 1.4301 – ASTM 304) is used, which stands out through its high stability, optimal hygienic properties and corrosion-resistance towards many (but not all!) chemical compounds (caution for example with chlorine compounds).

The chamber load for the appliance must be carefully checked for chemical compatibility with the materials mentioned. A material resistance table can be requested from the manufacturer.

## 2.5 Electrical equipment

- ▶ Operating voltage and current consumption: See nameplate
- ▶ Protection class I, i.e. operating insulation with PE conductor in accordance with EN 61010
- ▶ Protection type IP 20 acc. to EN 60 529
- ▶ Interference suppression acc. to EN 55011 class B
- ▶ Appliance fuse: Fusible link 250 V/15 A quick-blow
- ▶ The temperature controller is protected with a miniature fuse 100 mA (160 mA at 115 V)



## 2.6 Connections and interfaces

### 2.6.1 Electrical connection

This appliance is intended for operation on an electrical power system with a system impedance  $Z_{\max}$  of a maximum of 0.292 ohm at the point of transfer (service line). The operator must ensure that the appliance is operated only on an electrical power system that meets these requirements. If necessary, you can ask your local energy supply company what the system impedance is.

Observe the country-specific regulations when connecting (e.g. in Germany DIN VDE 0100 with residual current circuit breaker).

### 2.6.2 Communication interfaces

The interfaces are intended for appliances which meet the requirements of IEC 60950-1.

#### USB interface

The appliance is fitted by default with a USB interface in accordance with the USB specification. This way, you can

- ▶ transfer software stored on a USB storage medium to the appliance (see page 52).
- ▶ export protocol logs from the appliance to a USB storage medium (see page 54).
- ▶ transfer user ID data stored on a USB storage medium to the appliance (see page 55).

The USB interface is located on the lower right of the ControlCOCKPIT (Fig. 4).

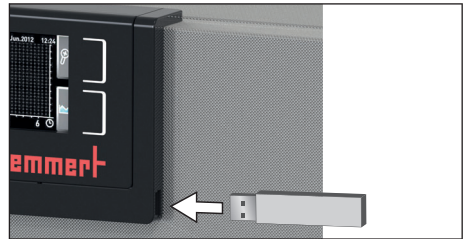


Fig. 4 USB interface

#### Ethernet interface

Via Ethernet interface, the appliance can be connected to a network, so that you can transfer programmes created with AtmoCONTROL software to the appliance and read out protocol logs. The Ethernet interface is located on the rear of the appliance (Fig. 5).

For identification purposes, each appliance connected must have its own unique IP address. Setting the IP address is described on page 44.

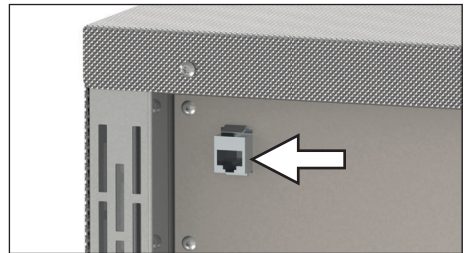


Fig. 5 Ethernet interface



You will find a description of how to transfer programmes via Ethernet in the enclosed AtmoCONTROL manual.

With an optional USB to Ethernet converter, the appliance can be directly connected to a computer / laptop (see Scope of delivery on page 16).

## 2.7 Designation (nameplate)

The nameplate (Fig. 6) provides information about the appliance model, manufacturer and technical data. It is attached to the front of the appliance, on the right side behind the door (see page 9).

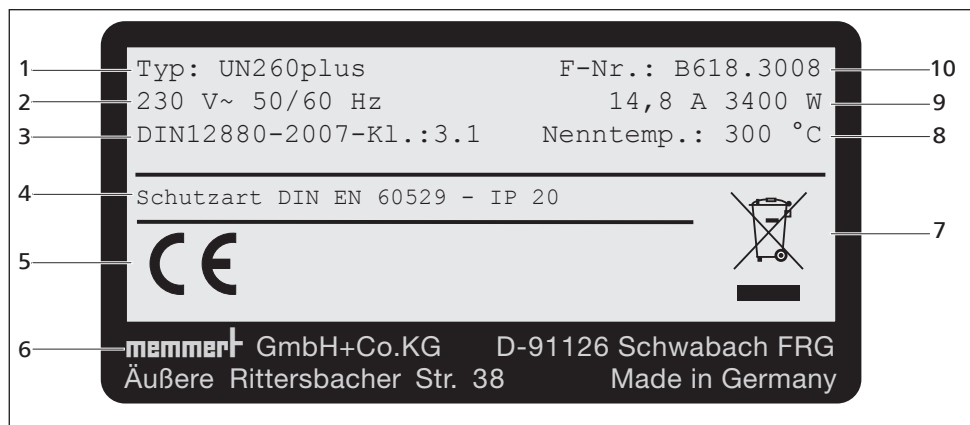


Fig. 6 Nameplate (example)

- |   |                   |    |                            |
|---|-------------------|----|----------------------------|
| 1 | Type designation  | 6  | Address of manufacturer    |
| 2 | Operating voltage | 7  | Disposal note              |
| 3 | Applied standard  | 8  | Temperature range          |
| 4 | Protection type   | 9  | Connection / power ratings |
| 5 | CE conformity     | 10 | Appliance number           |

## 2.8 Technical data

Appliance size		30	55	75	110	160	260	
Appliance width D <sup>1</sup> [mm]		585	585	585	745	745	824	
Appliance height E <sup>1</sup> [mm]		707	787	947	867	1107	1186	
Appliance depth F <sup>1</sup> (footprint) [mm]		434	514	514	584	584	684	
Depth of door lock [mm]		56						
Working chamber width A <sup>1</sup> [mm]		400	400	400	560	560	640	
Working chamber height B <sup>1</sup> [mm]		320	400	560	480	720	800	
Working chamber depth C <sup>1</sup> [mm]		250	330	330	400	400	500	
Chamber volume [litres]		32	53	74	108	161	256	
Weight [kg]		48	57	66	78	96	110	
Power [W]	I...	115 V, 50/60 Hz	800	900	900	900	900	900
		230 V, 50/60 Hz	1600	1000	1250	1400	1600	1700
	U...	230 V, 50/60 Hz	1600	2000	2500	2800	3200	3400
		115 V, 50/60 Hz	1600	1700	1800	1800	1800	1800
Current consumption [A]	I...	230 V, 50/60 Hz	7,0	4,3	5,4	6,1	7,0	7,4
		115 V, 50/60 Hz	7,0	7,8	7,8	7,8	7,8	7,8
	U...	230 V, 50/60 Hz	7,0	8,7	10,9	12,2	13,9	14,8
		115 V, 50/60 Hz	13,9	14,8	15,6	15,6	15,6	15,6
max. number of sliding shelves		3	4	6	5	8	9	
max. load per sliding shelf [kg]		20						
max. load per appliance [kg]		60	80	120	175	210	300	
Setting temperature range	IN/IF	+20 bis +80 °C <sup>2</sup>						
	UN/UF	+20 bis +300 °C <sup>3</sup>						
Adjustment precision	IN/IF	0,1 K						
	UN/UF	bis 99,9 °C: 0,1 K, ab 100 °C: 0,5 K						

<sup>1</sup> See Fig. 7 on page 15

<sup>2</sup> With the interior lighting on, the minimum temperature might not be reached.

<sup>3</sup> to +250 °C if your device is equipped with a glass door.

Appliance size		450	750	1060
Appliance width D <sup>1</sup> [mm]		1224	1224	1224
Appliance height E <sup>1</sup> [mm]		1247	1720	1720
Appliance depth F <sup>1</sup> (footprint) [mm]		784	784	1035
Depth of door lock [mm]		56		
Working chamber width A <sup>1</sup> [mm]		1040	1040	1040
Working chamber height B <sup>1</sup> [mm]		720	1200	1200
Working chamber depth C <sup>1</sup> [mm]		600	600	850
Chamber volume [litres]		449	749	1060
Weight [kg]		170	217	252

Appliance size		450	750	1060	
Power [W]	I...	115 V, 50/60 Hz	1500	1800	–
		230 V, 50/60 Hz	1800	2000	–
	U...	400 V, 50/60 Hz	5800	7000	7000
		3 x 230 V o. N., 50/60 Hz	5800	7000	7000
		3 x 208 V, 50/60 Hz	4800	5700	5700
Current consumption [A]	I...	230 V, 50/60 Hz	7,8	8,7	–
		115 V, 50/60 Hz	13,0	15,6	–
	U...	400 V, 50/60 Hz	3 x 8,4	3 x 10,1	3 x 10,1
		3 x 230 V o. N., 50/60 Hz	3 x 14,6	3 x 17,6	3 x 17,6
		3 x 208 V, 50/60 Hz	3 x 13,3	3 x 15,1	3 x 15,1
max. number of sliding shelves		8	14	14	
max. load per sliding shelf [kg]		30		60	
max. load per appliance [kg]		300			
Setting temperature range	IN/IF	+20 bis +80 °C <sup>2</sup>		–	
	UN/UF	+20 bis +300 °C <sup>2 3</sup>		–	
Adjustment precision	IN/IF	0,1 K		–	
	UN/UF	bis 99,9 °C: 0,1 K, ab 100 °C: 0,5 K			

<sup>1</sup> See Fig. 7 on page 15

<sup>2</sup> With the interior lighting on, the minimum temperature might not be reached.

<sup>3</sup> to +250 °C if your device is equipped with a glass door.

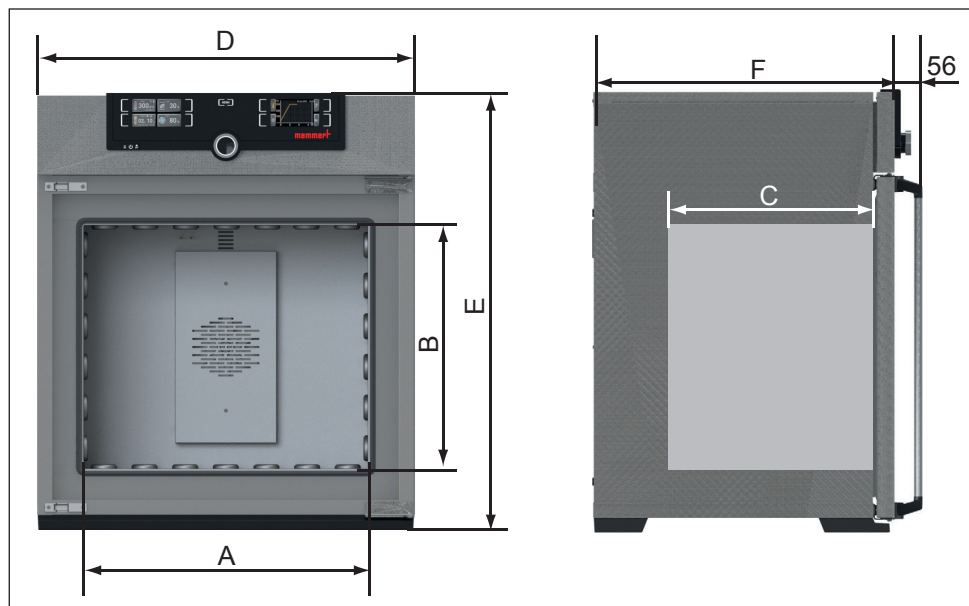


Fig. 7 Dimensions (see table on page 14)

## 2.9 Declaration of conformity

You can download the EC declaration of conformity of the appliance online:

English: <http://www.memmert.com/en/service/downloads/ce-statement/>

German: <http://www.memmert.com/de/service/downloads/eg-konformitaetserklaerung/>

## 2.10 Ambient conditions

- ▶ The appliance may only be used in enclosed rooms and under the following ambient conditions:

Ambient temperature	+5 °C to +40 °C
Humidity rh	max. 80 %, non-condensing
Overvoltage category	II
Pollution degree	2
Altitude of installation	max. 2,000 m above sea level

- ▶ The appliance may not be used in areas where there is a risk of explosion. The ambient air must not contain any explosive dusts, gases, vapours or gas-air mixtures. The appliance is not explosion-proof.
- ▶ Heavy dust production or aggressive vapours in the vicinity of the appliance could lead to sedimentation in the interior and, as a consequence, could result in short circuits or damage to electrical parts. For this reason, sufficient measures to prevent large clouds of dust or aggressive vapours from developing should be taken.

## 2.11 Scope of delivery

- ▶ Power cable
- ▶ Tilt protection
- ▶ One or two sliding steel grids (load capacity 30 kg each)
- ▶ USB storage medium with software and AtmoCONTROL manual
- ▶ The operating instructions at hand
- ▶ Calibration certificate

## 2.12 Optional accessories

- ▶ USB to Ethernet converter (Fig. 8). Makes it possible to connect the appliance's network interface (see page 12) to the USB port of a computer / laptop.
- ▶ Reinforced, sliding steel grids with a load capacity of 60 kg each (for appliance size 110 and larger)

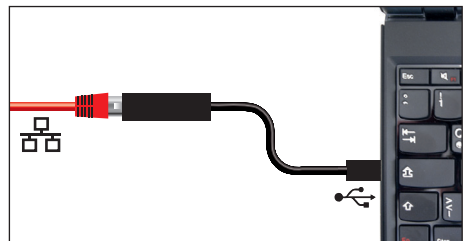


Fig. 8 Converter USB to Ethernet

### 3. Delivery, transport and setting up

#### 3.1 For your Safety



**Warning!**

Because of the heavy weight of the appliance, you could injure yourself if you try to lift it. To carry appliances of the sizes 30 and 55, at least two persons, for appliances of the sizes 75, 110, 160 and 260, four people are needed. Appliances larger than that may not be carried but must be transported with a manual pallet jack or forklift truck.

30	55	75	110	160	260	450	750



**Warning!**

You may get your hands or feet squashed when transporting and installing the appliance. Wear protective gloves and safety boots. When grasping the bottom of the appliance, grasp it only on the sides:



**Warning!**

The appliance could fall over and seriously injure you. Never tilt the appliance and transport it in upright position and without load only (except for standard accessories such as steel grids or shelves). Appliances with castors always have to be moved by two people.

## 3.2 Delivery

The appliance is packed in cardboard and is delivered on a wooden palette.

## 3.3 Transport

The appliance can be transported in three ways:

- ▶ With a forklift truck; move the forks of the truck entirely under the pallet
- ▶ On a manual pallet jack
- ▶ On its own castors, in case of the corresponding configuration, for which the catch on the (front) castors must be released

## 3.4 Unpacking

**i** To avoid damage, do not unpack the appliance until you reach the installation site.

Remove the cardboard packaging by pulling it upwards or carefully cutting along an edge.

### 3.4.1 Checking for completeness and transport damage

- ▶ Check the delivery note to ensure that the delivery is complete.
- ▶ Check the appliance for damage.

If you notice deviations from the delivery note, damage or irregularities, do not put the appliance into operation but inform the haulage company and the manufacturer.

### 3.4.2 Removing the transport protection

Remove the transport protection. It is located between the door hinge, door and frame and has to be removed after opening the door.

### 3.4.3 Disposing of packaging material

Dispose of the packaging material (cardboard, wood, foil) in accordance with the applicable disposal regulations for the respective material in your country.

## 3.5 Storage after delivery

If the appliance is first to be stored after delivery: Read the storage conditions from page 57.



## 3.6 Setting up



### Warning!

Due to its centre of gravity, the appliance can fall over to the front and injure you or other people. Always attach the appliance to a wall with the tilt protection (see page 21). If this cannot be done due to space problems, do not operate the appliance and do not open the door. Contact the Memmert service team (see page 2).

### 3.6.1 Prerequisites

The installation site must be flat and horizontal and must be able to reliably bear the weight of the appliance (see "Technical data" on page 14). Do not place the appliance on a flammable surface.

Depending on the model (see nameplate), a 230 V, 115 V or 400 V power connection must be available at the installation site.

The distance between the wall and the rear of the appliance must be at least 15 cm. The clearance from the ceiling must not be less than 20 cm and the side clearance from walls or nearby appliances must not be less than 5 cm (Fig. 9). Sufficient air circulation in the vicinity of the appliance must be guaranteed at all times.

For appliances with castors, these need to be positioned in forward direction at all times.

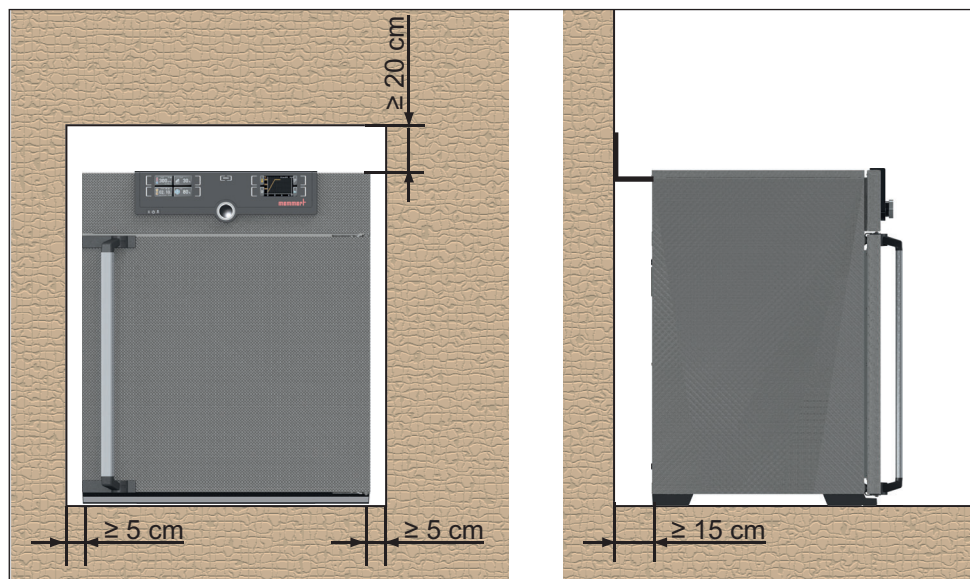


Fig. 9 Minimum clearance from walls and ceiling



3.6.2 Installation options

Setting up	Comments	Suitable for appliance size ...							
		30	55	75	110	160	260	450	750
<p>Floor</p>		✓	✓	✓	✓	✓	✓	✓	✓
<p>Table</p>	Check the load capacity first	✓	✓	✓	✓	✗	✗	✗	✗
<p>Stacked</p>	two appliances maximum; mounting material (feet) provided	✓	✓	✓	✓	✗	✗	✗	✗
<p>Wall mounting</p>	Separately packaged fastening material is included in the scope of delivery. Observe the assembly instructions provided.	✓	✓	✓	✓	✓	✗	✗	✗
<p>Base</p>	with/without castors	✓	✓	✓	✓	✓	✓	✓	✗
<p>Castor frame</p>		✓	✓	✓	✓	✓	✓	✗	✗
<p>Height adjustable feet</p>		✓	✓	✓	✓	✓	✓	✓	✓

### 3.6.3 Tilt protection

Attach the appliance to a wall with the tilt protection. The tilt protection is included in the delivery.

1. As illustrated, fasten the tilt protection to the rear side of the appliance.
2. Bend the tilt protection upwards by  $90^\circ$  in the desired distance to the wall (consider the minimum distance to the wall, see Fig. 9).
3. Drill a hole, insert a plug and screw the tilt protection to a suitable wall.

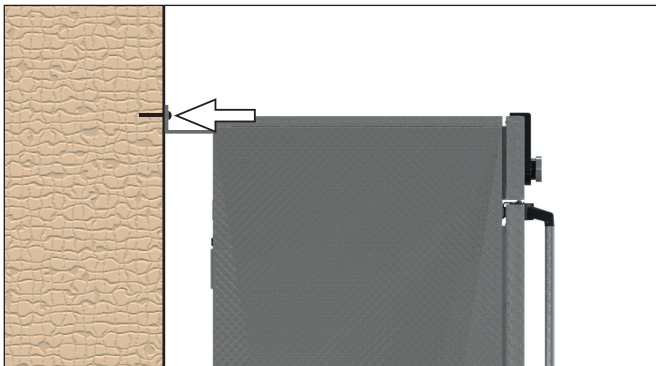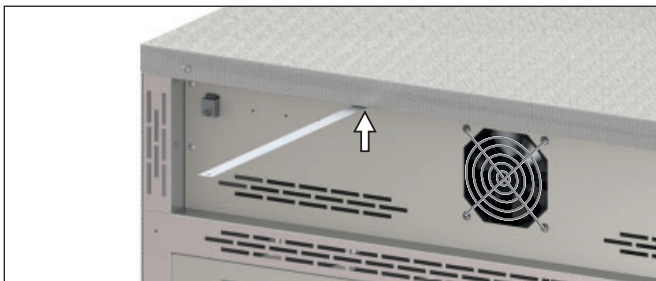

### 3.6.4 Adjusting the doors

For appliances it is possible to adjust doors that warp due to the floor conditions. In order to do so, every door has two adjuster screws at the top and at the bottom (Fig. 10).

**1** First, adjust the door at the top and then, if further adjustment is necessary, at the bottom as well.

1. Open the door.
2. Undo the screws.
3. Adjust the door.
4. Tighten the screws again.
5. Check door alignment.
6. If necessary, readjust.

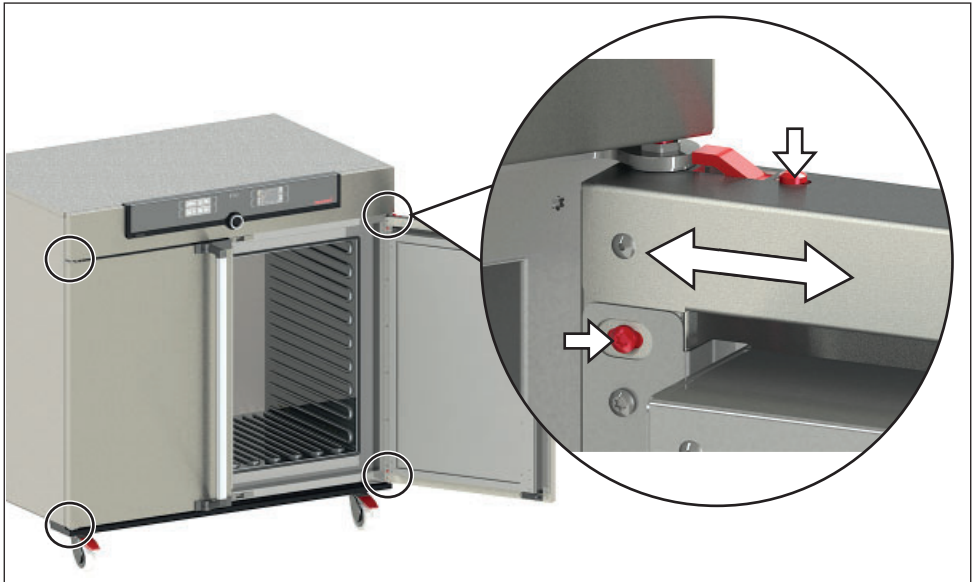


Fig. 10 Door adjustment screws

## 4. Putting into operation

### Caution:

The first time the appliance is operated, it must not be left unattended until it has reached the steady state.

### 4.1 Connecting the appliance

#### ⚠ WARNING



**Condensation might cause a short circuit. After transport or storage at high humidity conditions, the appliance shall be stored unpacked at normal conditions for at least 24 hours. During this period of time the appliance shall not be connected to the power supply.**

### Caution:

Observe the country-specific regulations when making connections (e.g. in Germany DIN VDE 0100 with residual current circuit breaker). Observe the connection and power ratings (see nameplate and Technical data on page 14). Make sure to establish a safe PE conductor connection.

Lay the power cable so that

- ▶ it is always accessible and within reach so it can be disconnected quickly in the event of failure or emergencies;
- ▶ no one can trip over it;
- ▶ it does not come into contact with any hot parts.

#### 230/115-V appliances:

Plug the provided power cable into the rear of the appliance and connect it to a CEE 7/4 socket (Fig. 11).

#### 400 V appliances:

The power cable is permanently installed. Connect the plug to a 400 V CEE coupling (Fig. 12).

#### 3 x 208 V appliances:

The power cable is permanently installed. Connect the plug to a 3 x 208 V / 20 A coupling (NEMA L15-20R) (Fig. 13).

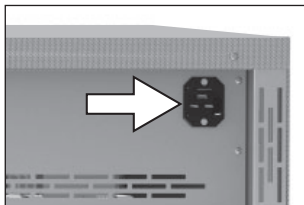


Fig. 11 Power connection  
230/115 V




Fig. 12  
400 V CEE connection



Fig. 13 NEMA L15-20R  
twist lock connection

## 4.2 Switching on

Switch on the appliance by pressing the On/Off switch on the front of the appliance (Fig. 14).

The starting process is shown by three animated white dots . If the dots have another colour, an error has occurred (see page 40).

**i** After the first start-up, the appliance display is set to English by default. You can change the language as described from page 43. However, to get a basic overview of operating the appliance, you should read the following chapter first.

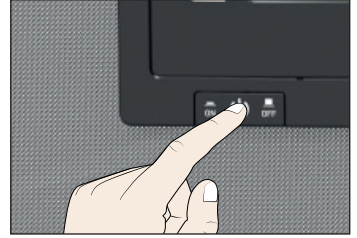


Fig. 14 Switch on appliance

## 5. Operation and control

### 5.1 Operating personnel

The appliance may only be operated by persons who are of legal age and have been instructed accordingly. Personnel who are to be trained, instructed or who are undergoing general training may only work with the appliance under the continuous supervision of an experienced person.

### 5.2 Opening the door

- ▶ To open the door, pull the door handle to the side (to the left or to the right, depending on the door variation, see Fig. 15, A) and open the door completely.
- ▶ To close the appliance, push the door closed and the door handle to the side (B).

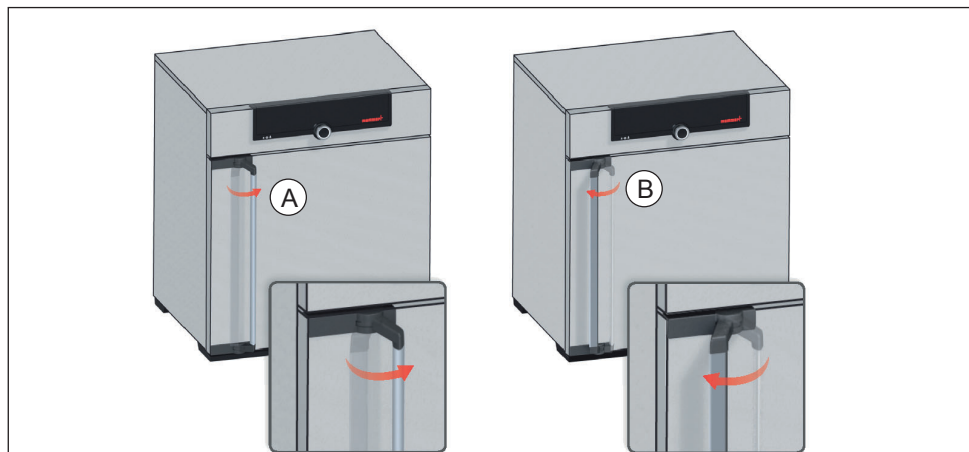


Fig. 15 Opening and closing the door



#### Warning!

If the door is open while the appliance is in operation, the appliance may overheat and pose a fire hazard. Do not leave the door open during operation.



#### Warning!

In case of appliances of a certain size, you can get accidentally locked in, which is life-threatening. Do not climb into the appliance!



## 5.3 Loading the appliance



### Warning!

When loading the appliance with an unsuitable load, poisonous or explosive vapours or gases may be produced. This could cause the appliance to explode, and persons could be severely injured or poisoned. The appliance may only be loaded with materials which do not form any toxic or explosive vapours when heated up and cannot ignite (see also page 10). If there is any doubt as to the composition of materials, they must not be loaded into the appliance.



### i Caution:

Check the chamber load for chemical compatibility with the materials of the appliance (see page 11).

### i Caution:

If the chamber load is wet or very humid, water may accumulate on the floor and damage the heater. If wet, use a drip tray on the bottom of the tube.

Insert the sliding steel grids or sliding shelves. The maximum number of grids / shelves and the load capacity are specified in the technical data overview from page 14.

The chamber must not be loaded too tightly, so that proper air circulation in the working chamber is guaranteed. Do not place any chamber load on the floor, touching the side walls or right below the ceiling of the working chamber (Fig. 16, see also the "correct loading" sticker on the appliance).

In case of improper loading (chamber loaded too tightly), the set temperature may be exceeded or it may take longer until it is reached.

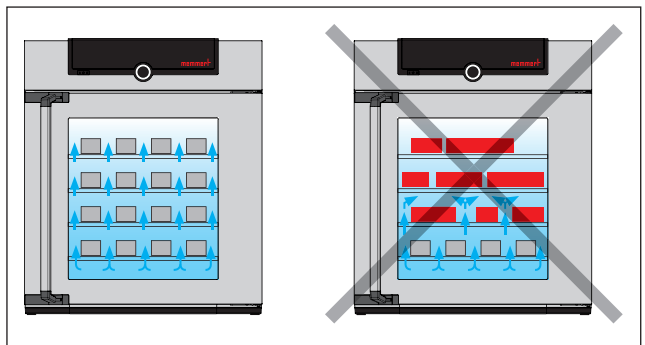


Fig. 16 Correct placement of the chamber load

i To achieve the correct heating capacity, the type of slide-in unit used – Grid or Shelf – must be set in the menu under SETUP (see page 44).

## 5.4 Operating the appliance

### 5.4.1 ControlCOCKPIT

In manual operation, the desired parameters are entered at the ControlCOCKPIT on the front of the appliance (Fig. 17 and Fig. 18). You can also make basic settings here (menu). Additionally, warning messages are displayed, e.g. if the temperature is exceeded. In programme mode, the parameters defined, the programme description, the programme segment currently active and programme duration remaining are displayed (for a more detailed description, see page 31).

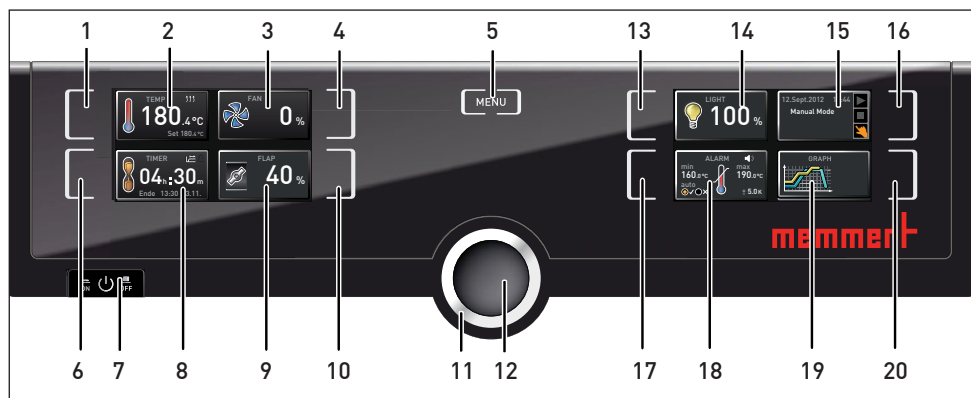


Fig. 17 ControlCOCKPIT for UF.../IF... appliances in operating mode (width may differ depending on appliance size)

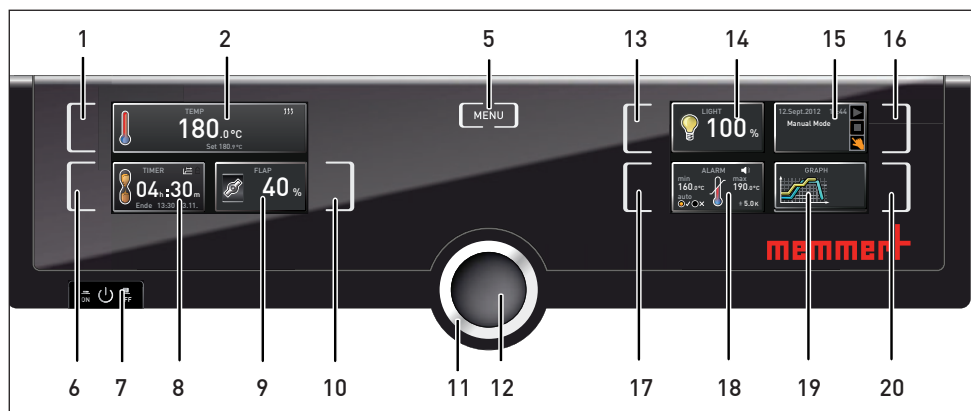



Fig. 18 ControlCOCKPIT for UN.../IN... appliances in operating mode (width may differ depending on appliance size)

- |  |   |
|--|---|
| <ol style="list-style-type: none"> <li>1 Activation key for temperature setpoint adjustment</li> <li>2 Setpoint and actual temperature display</li> <li>3 Fan speed display</li> <li>4 Activation key for fan speed setting</li> <li>5 Switch to menu mode (see page 42)</li> <li>6 Activation key digital backwards counter with target time setting, adjustable from 1 minute to 99 days</li> <li>7 On/Off switch</li> <li>8 Display digital backwards counter with target time setting, adjustable from 1 minute to 99 days</li> <li>9 Air flap position display</li> </ol> | <ol style="list-style-type: none"> <li>10 Activation key for air flap position adjustment</li> <li>11 Turn control for setpoint adjustment</li> <li>12 Confirmation key (accepts setting made with the turn control)</li> <li>13 Activation key for interior lighting (additional option)</li> <li>14 Interior lighting display (additional option)</li> <li>15 Appliance state and programme display</li> <li>16 Activation key for the appliance state</li> <li>17 Activation key for temperature monitoring</li> <li>18 Temperature monitoring display</li> <li>19 Graphic representation</li> <li>20 Activation key for graphic representation</li> </ol> |
|--|---|


### 5.4.2 Basic operation

In general, all settings are made according to the following pattern:


1. Activate the desired parameter (e.g. temperature). To do so, press the corresponding activation key on the left or right or the respective display. The activated display is lined in colour, the other displays are dimmed. The set value is highlighted in colour.

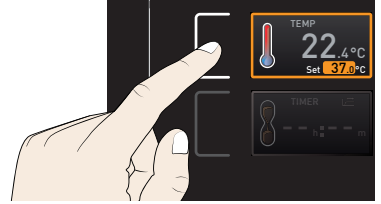

2. By turning the turn control to the left or right, adjust the set value (e.g. to 180.0 °C).

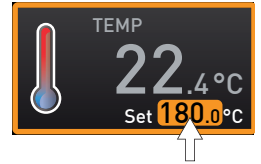


3. Save the set value by pressing the confirmation key. The display returns to normal and the appliance begins adjusting to the defined set value.

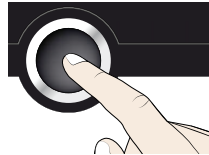

Additional parameters (air flap position etc.) can be set accordingly.

- 1 If no new values are entered or confirmed for approx. 30 seconds, the appliance automatically returns to the main menu and restores the former values.

If you want to cancel the setting procedure, press the activation key on the left or right of the display that you want to exit. The appliance restores the former values. Only the settings that you have confirmed by pressing the confirmation key before cancelling the setting procedure are accepted.

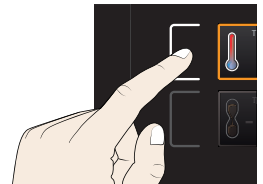

### 5.4.3 Operating modes



The appliance can be operated in different modes:

- ▶ **Manual mode:** The appliance runs in permanent operation at the values set on the ControlCOCKPIT. Operation in this mode is described in chapter 5.4.4.
- ▶ **Operation with digital backwards counter with target time setting, adjustable from 1 minute to 99 days (Timer):** The appliance runs at the values set until the timer has elapsed. Operation in this mode is described in chapter 5.4.5.
- ▶ **Programme mode:** The appliance automatically runs programme sequences which have been defined using AtmoCONTROL software at a computer / laptop and then transferred to the appliance from a USB stick or via Ethernet. Operation in this mode is described in chapter 5.4.6.
- ▶ via remote control


**i** The status display shows you which operating mode or operating state the appliance is currently in. The current operating state is highlighted in colour and indicated by the text display:

- ▶ Appliance is in programme mode
- Programme is stopped
- ✎ Appliance is in manual operating mode

The example on the right shows the appliance in manual mode, identified by the coloured hand symbol.

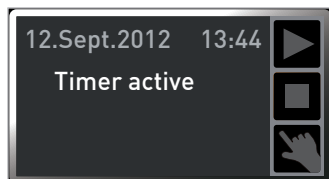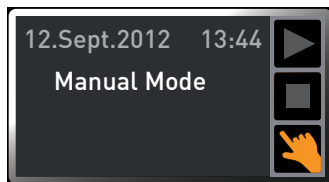

▶ When the appliance is in timer mode, Timer active is displayed:

▶ When the appliance is in remote control mode, the  symbol appears in the temperature display:

### 5.4.4 Manual mode

In this operating mode, the appliance runs in permanent operation at the values set on the ControlCOCKPIT.

#### Adjustment options



As described in chapter 5.4.2, you can set the following parameters after pressing the corresponding activation key (in any sequence):

#### Temperature

Adjustment range: model dependent (see nameplate and technical date on page 14)


**i** Heating operation is indicated by the  symbol.

**i** You can select °C or °F as the temperature unit displayed (see page 45).

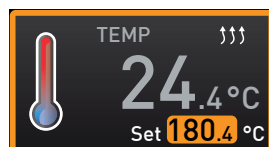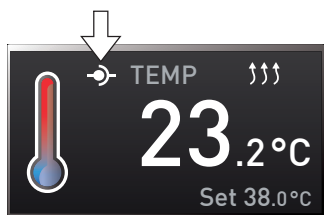

#### Air flap position

Adjustment range: 0 % (closed, recirculating operation) to 100 % (completely opened, fresh air operation) in steps of 10 %

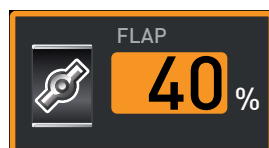

#### Fan speed


(only for UF.../IF... appliances)

Adjustment range: 0 to 100 % in steps of 10%



#### Interior lighting (additional option)


Adjustment range: 0 %, 100 %

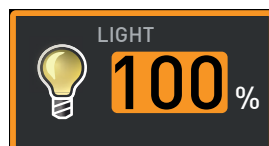

### 5.4.5 Operation with digital backwards counter with target time setting, adjustable from 1 minute to 99 days (Timer)

In timer operation, you can adjust the time the appliance runs at the set value. The appliance has to be in manual operating mode for this.

1. Press the activation key to the left of the timer display. The timer display is activated.

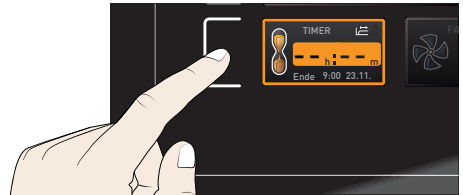


2. Turn the turn control until the desired duration is displayed – in this example 4 hours 30 minutes. The approximate end time is shown beneath, in a smaller font.




- i** Up to a duration of 23 hours 59 minutes, the time is displayed in hh:mm (hours:minutes) format. For 24 hours and more, the format dd:hh (days:hours) is used. The maximum duration adjustable is 99 days 00 hours.


3. Press the confirmation key to confirm.

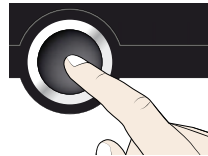


The display now shows the remaining time in a large font and the approximate end time in a smaller font beneath. The status display shows „Timer active“.

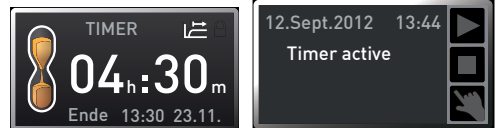


4. Now, as described under 5.4.2, set the individual values for temperature, air flap position etc. which you want the appliance to operate at. The set values can be changed while the timer elapses. The changes are effective immediately.

- i** In Setup, you can choose if the timer should run setpoint-dependent or not. This determines whether the timer should not start until a tolerance band around the set temperature is reached or whether it should start right after activation (see page 46). If the timer runs setpoint-dependent, this is indicated by the  symbol in the timer display.

When the timer has elapsed, the display shows 00h:00m. All functions (heating etc.) are switched off. If a fan had been active, it will keep on running for a short safety period. In addition, an acoustic alarm sounds, which can be turned off by pressing the confirmation key.

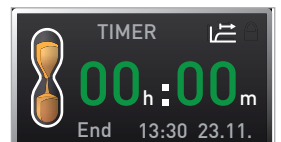

To deactivate the timer, open the timer display by pressing the activation key again and then turning the turn control to reduce the timer setting until --:-- is displayed. Confirm with the confirmation key.

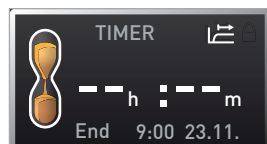

### 5.4.6 Programme mode

In this operating mode, programmes saved in the appliance can be started with different combinations of individual parameters (temperature, air flap position, fan speed, working chamber lighting) at staggered intervals, which the appliance then automatically processes in sequence. These programmes are not created directly at the appliance but externally at a computer / laptop and using AtmoCONTROL software. Transfer to the appliance is possible using the provided USB storage medium or via Ethernet.




A description of how to create and save programmes can be found in the separate AtmoCONTROL software manual.

#### Starting a programme

1. Press the activation key to the right of the status display. The current operating mode is highlighted automatically, in this example Manual Mode (👉).




2. Turn the turn control until the ▶ start symbol is highlighted. The current programme is displayed, in this example Test 012.

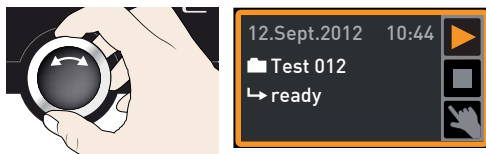

**i** Only the programme currently selected in the menu and shown in the display can be used. If you want to process another programme, you need to activate it in the menu first (description from page 48).

3. To start the programme, press the confirmation key. The programme starts. The display shows:


- ▶ the programme description (in this example Test 012)
- ▶ the programme segment description, in this example Ramp 1
- ▶ the current run (in case of loops)



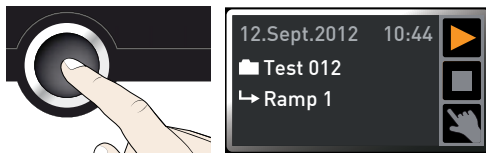



**i** You cannot change any parameters (e.g. the temperature) at the appliance while a programme is running. However, the displays ALARM and GRAPH can still be used.

### Canceling a programme

You can cancel an active programme at any time.


1. Press the activation key to the right of the status display. The status display is automatically highlighted.
2. Turn the turn control until the stop symbol  is highlighted.
3. Press the confirmation key to confirm. The programme is cancelled.

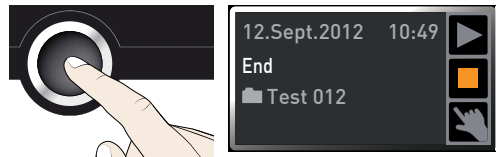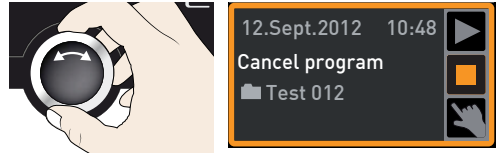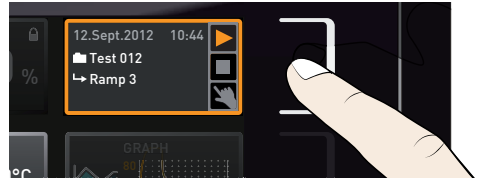


 A cancelled programme cannot be resumed at the point it was cancelled. It must be restarted from the beginning.


### End of programme

The display shows End when the programme is finished.

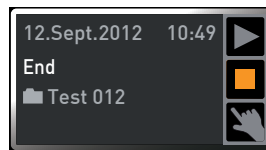

You can now

- ▶ restart the programme as described
- ▶ select another programme for processing in menu mode (see page 52) and run it as described.
- ▶ return to manual mode. To do so, reactivate it by pressing the activation key next to the status display, then turn the turn control until the hand symbol  is highlighted in colour and press the confirmation key.

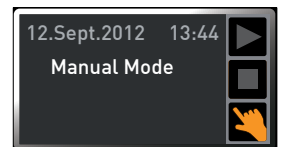




## 5.5 Temperature monitoring

The appliance is equipped with a multiple overtemperature protection (mechanical/electronic) in accordance with DIN 12 880. This serves to avoid damage to the chamber load and/or appliance in case of a malfunction:

- ▶ electronic temperature monitoring (TWW/TWB)
- ▶ automatic temperature monitor (ASF)
- ▶ mechanical temperature limiter (TB)

The monitoring temperature of the electronic temperature monitoring is measured via a separate Pt100 temperature sensor in the working chamber. Temperature monitoring settings are made via the ALARM display. The settings made apply to all operating modes.

If temperature monitoring has been triggered, this is indicated by the temperature display: the actual temperature is highlighted in red and a warning symbol  is shown (Fig. 19). The type of temperature monitoring triggered (TWW in this example) is shown beneath the temperature. If the acoustic alarm has been activated in the menu mode (Sound, see page 53, indicated by the speaker symbol in the alarm display), the alarm is additionally signalled by an intermittent acoustic signal, which can be turned off by pressing the confirmation key. Information on what to do in this case are provided in the chapter Malfunctions, warning and error messages from page 38.

Before reading how to adjust temperature monitoring (from page 35), please read the description of the individual monitoring functions here.

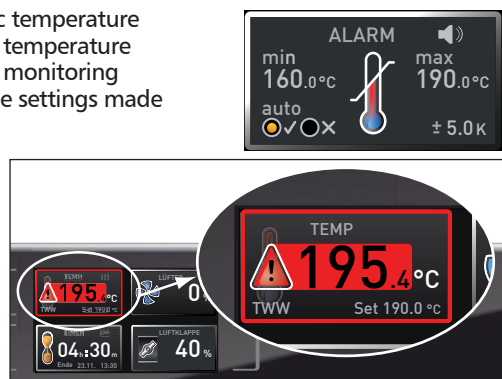


Fig. 19  
Temperature monitoring triggered

5.5.1 Electronic temperature monitoring (TWW)

The manually set monitoring temperature min and max of the electronic overtemperature control is monitored by an adjustable over/undertemperature controller (TWW) protection class 3.1 acc. to DIN 12 880 (or over/undertemperature controller (TWW) protection class 3.1 for UIS appliances). If the manually set monitoring temperature max is exceeded, the TWW takes over temperature control and begins to regulate the monitoring temperature (Fig. 20).

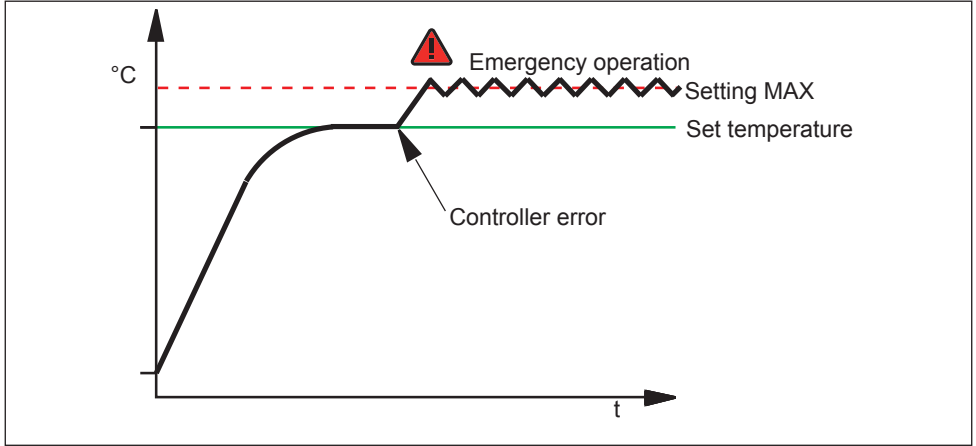


Fig. 20 Schematic diagram of how the TWW temperature monitoring works

5.5.2 Electronic temperature limiter (TWB) protection class 2 acc. to DIN 12 880

If the manually set monitoring temperature max is exceeded, the TWB switches off heating permanently (Fig. 21) and can be reset by pressing the confirmation key.

**i** In programme mode, the current programme is resumed for TWB alarms of up to 15 minutes. If the alarm is active for more than 15 minutes, the programme is cancelled.

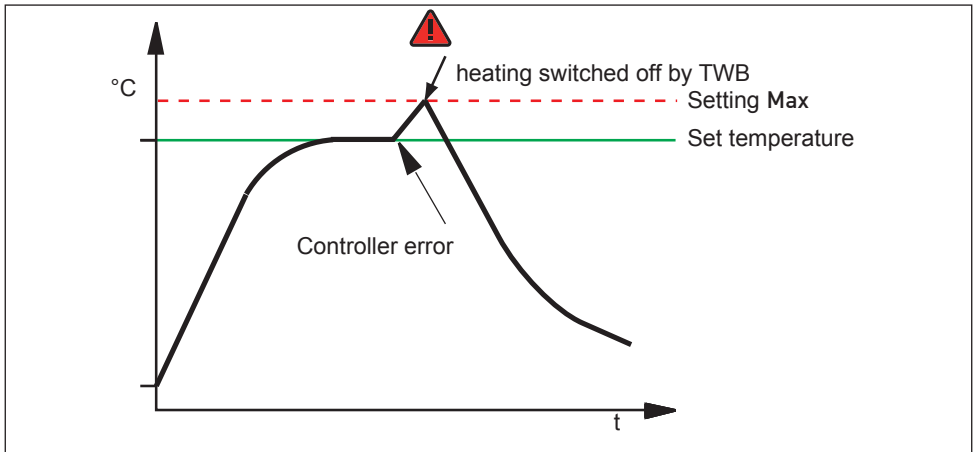


Fig. 21 Schematic diagram of how the TWB temperature monitoring works

### 5.5.3 Automatic temperature monitor (ASF)

ASF is a monitoring device that automatically follows the set temperature setpoint within an adjustable tolerance band (Fig. 22).

The ASF – if switched on – is automatically activated as soon as the actual temperature value reaches 50 % of the set tolerance band of the setpoint (in the example: 180 °C – 1.5 K) for the first time (section A).

When the temperature violates the set tolerance band around the setpoint (in the example in Fig. 22:

180 °C ± 3 K) – e.g. if the door is opened during operation (section B of illustration) – the alarm is set off. The ASF alarm is automatically triggered as soon as 50 % of the set tolerance band of the setpoint (in the example: 180 °C ± 1.5 K) are reached again (section C).

If the temperature setpoint is altered, the ASF is automatically disabled temporarily (in this example: The setpoint is changed from 180 °C to 173 °C, section D) until the tolerance range of the new temperature setpoint is reached again (section E).

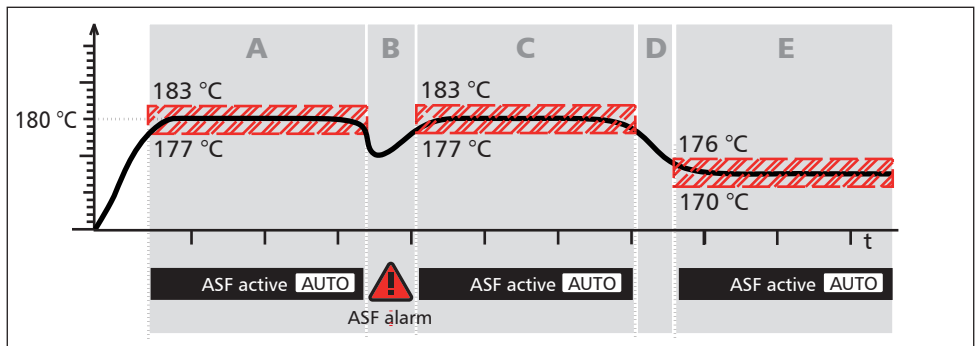


Fig. 22 Schematic diagram of how the ASF temperature monitoring works


### 5.5.4 Mechanical temperature monitoring: Temperature limiter (TB)

The appliance is equipped with a mechanical temperature limiter (TB) of protection class 1 in accordance with DIN 12 880.


If the electronic monitoring unit should fail during operation and the factory-set maximum temperature is exceeded by approx. 20 °C, the temperature limiter, as the final protective measure, switches off the heating permanently.

### 5.5.5 Adjusting temperature monitoring



1. Press the activation key to the left of the ALARM display. The min setting (under-temperature protection) is automatically activated.

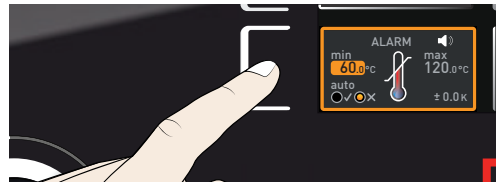

- By turning the turn control, adjust the desired lower alarm limit value, in the example on the right 160 °C.
- If no undertemperature protection limit is required, set the lowest temperature.

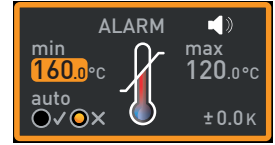


- Press the confirmation key to confirm. The max display (overtemperature protection) is activated.

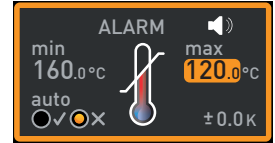
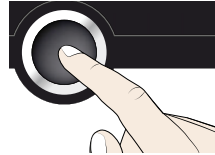


- By turning the turn control, adjust the desired upper alarm limit value, in the example on the right 190 °C.


- The monitoring temperature must be set sufficiently high above the maximum set temperature. We recommend 5 to 10 K.

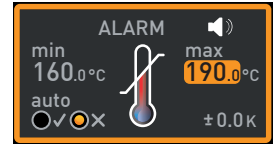



- Accept the upper alarm limit value by pressing the confirmation key. The setting of the automatic temperature monitor (ASF) is automatically activated (auto).

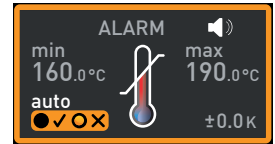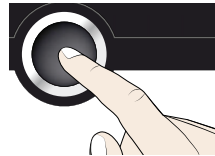

- With the turn control, select ON (✓) or OFF (✗).

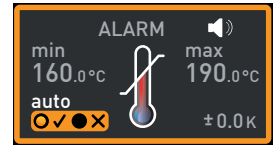


- Press the confirmation key to confirm. The ASF tolerance band setting is activated.

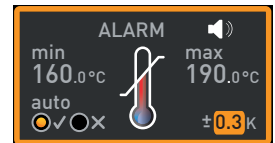
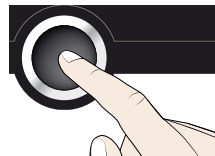


- With the turn control, adjust the desired tolerance band, e.g. 5.0 K.

- We recommend a tolerance band of 5 to 10 K (1 to 3 K for incubators IN/IF).

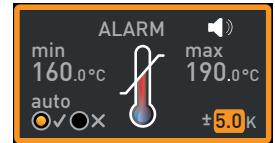

- Press the confirmation key to confirm. Temperature monitoring is now active.

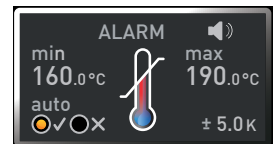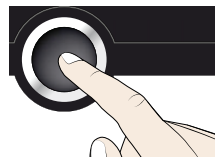


In the menu, you can set:

- ▶ which type of protection (TWW or TWB) should be active (see page 45)
- ▶ if an acoustic signal should be triggered on alarm (see page 53)


## 5.6 Graph

The GRAPH display provides an overview of the chronological sequence of the set values and the actual values as a curve.




1. Press the activation key to the right of the GRAPH display. The display is enlarged and the temperature profile shown.

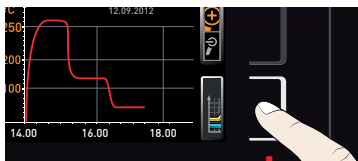

- ▶ To change the time frame to be displayed: Press the activation key next to the <> arrow symbols. The time frame to be displayed can now be changed by turning the turn control.

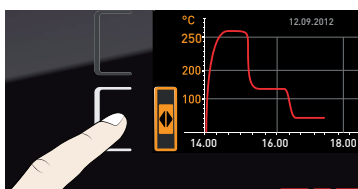

- ▶ To zoom the graph in or out: Press the activation key next to the magnifying glass symbol. Select whether you want to zoom in or out (+/-) with the turn control and confirm your selection by pressing the confirmation key.

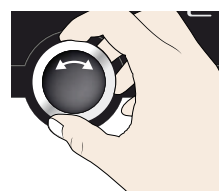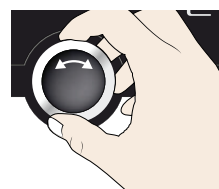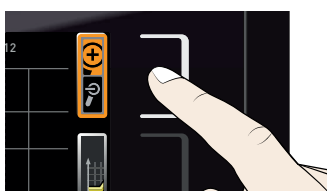

To close the graphical representation, again press the activation key which you have used to activate it.

## 5.7 Ending operation



### Warning!

Depending on the operation performed, the surfaces in the working chamber and the chamber load may still be very hot after the appliance is switched off. Touching these surfaces can cause burns. Wear heat-resistant protective gloves or wait until the appliance cools down.

1. Switch off active appliance functions (turn back the set values).
2. Remove the chamber load.
3. Switch off the appliance (Fig. 23).



Fig. 23 Switch off appliance

## 6. Malfunctions, warning and error messages




### Warning!


After removing covers, live parts may be exposed. You may receive an electric shock if you touch these parts. Malfunctions requiring work inside the appliance may only be rectified by electricians. Observe the separate service manual for this.

Do not try to rectify appliance errors yourself but contact the MEMMERT customer service department (see page 2) or an authorised service point.



In case of enquiries, please always specify the model and appliance number from the nameplate (see page 13).



### 6.1 Warning messages of the monitoring function

**1** If the acoustic alarm has been activated in the Signals menu (see page 53), which is indicated by the speaker symbol  in the alarm display, the alarm is additionally signalled by an intermittent acoustic signal. If the confirmation key is pressed, the acoustic alarm can be temporarily switched off until the next alarm event occurs.

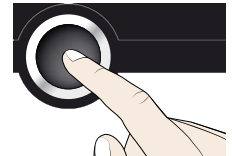





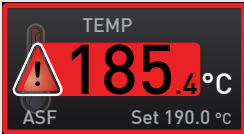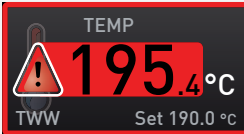

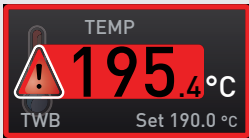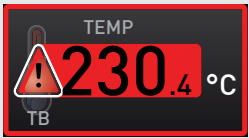
#### 6.1.1 Temperature monitoring

Description	Cause	Action	See
Temperature alarm and "ASF" are displayed 	Automatic temperature monitor (ASF) triggered	Check if the door is closed. Close the door. Extend the ASF tolerance band If the alarm continues: Contact customer service	page 36 page 2
Temperature alarm and "TWW" are displayed 	The adjustable undertemperature / overtemperature controller (TWW) has assumed heating control.	Increase the difference between the monitoring and setpoint temperature – by either increasing the max value of the temperature monitoring or decreasing the setpoint temperature. If the alarm continues: Contact customer service	page 35 page 2

Description	Cause	Action	See
Temperature alarm and "TWB" are displayed 	The electronic temperature limiter (TWB) permanently switched off heating.	Deactivate the alarm by pressing the confirmation key.  Increase the difference between the monitoring and setpoint temperature – by either increasing the max value of the temperature monitoring or decreasing the setpoint temperature.  If the alarm continues: Contact customer service	page 35   page 2
Temperature alarm and "TB" are displayed 	The mechanical temperature limiter (TB) permanently switched off heating.	Switch off the appliance and leave to cool down. Contact customer service and have the error rectified (e.g. by replacing the temperature sensor).	page 2
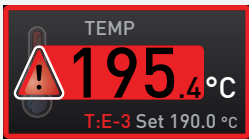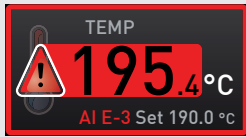
## 6.2 Malfunctions, operating problems and appliance errors

Error description	Cause of error	Rectifying errors	See
Displays are dark	External power supply was interrupted	Check the power supply	page 23
	Miniature fuse, appliance fuse or power module faulty	Contact customer service	page 2
Displays cannot be activated	Appliance locked by user ID	Unlock with user ID	page 55
	The appliance is in programme, timer or remote control mode (mode "Write" or "Write + Alarm")	Wait until the end of the programme or timer mode or switch off the remote control	
Displays suddenly look different	Appliance is in "wrong" mode	Change to operating or menu mode by pressing the MENU key	

Error description	Cause of error	Rectifying errors	See
Display T:E-3 in the temperature display 	Temperature operating sensor defective. The monitoring sensor takes over the measurement function.	<ul style="list-style-type: none"> <li>▶ The appliance can be temporarily operated</li> <li>▶ Contact customer service as soon as possible</li> </ul>	page 2
Error message AI E-3 in the temperature display 	Temperature monitoring sensor defective. The operating sensor takes over the measurement function.	<ul style="list-style-type: none"> <li>▶ The appliance can temporarily be kept in service</li> <li>▶ Contact customer service as soon as possible</li> </ul>	page 2
Error message E-3 in the temperature display 	Operating and monitoring sensor defective	<ul style="list-style-type: none"> <li>▶ Switch off appliance.</li> <li>▶ Remove the chamber load</li> <li>▶ Contact customer service</li> </ul>	page 2
When switching on the appliance, the start animation is displayed in another colour than white 	<ul style="list-style-type: none"> <li>▶ Cyan : Not enough storage space on the SD card</li> <li>▶ Red : The system files could not be loaded</li> <li>▶ Orange : The fonts and images could not be loaded</li> </ul>	Contact customer service  Contact customer service  Contact customer servi	page 2  page 2  page 2



### 6.3 Power failure

In case of a power failure, the appliance operates as follows:

#### In manual mode

After power supply has been restored, operation is continued with the parameters set. The time and duration of the power failure are documented in the log memory.

#### In timer or programme mode

In case of an interruption of the power supply of less than 60 minutes, the current programme is continued from the point at which it was interrupted. For interruptions of the power supply longer than this, all appliance functions (heating, fan etc.) are switched off and the air flap is opened.

#### In remote control mode:

The previous values are restored. If a programme has been initiated via remote control, it is continued.

## 7. Menu mode


In menu mode, you can make basic settings, load programmes and export protocols, as well as adjust appliance parameters.

### Caution:

**1** Before changing menu settings, read the description of the respective functions on the following pages to avoid possible damage to the appliance and/or chamber load.

To enter menu mode, press the MENU key.

**1** To exit the menu mode at any time, press the MENU key again. The appliance then returns to manual mode. Only changes accepted by pressing the confirmation key are saved.

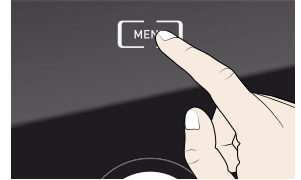

### 7.1 Overview

Press the MENU key to change between the displays in menu mode:

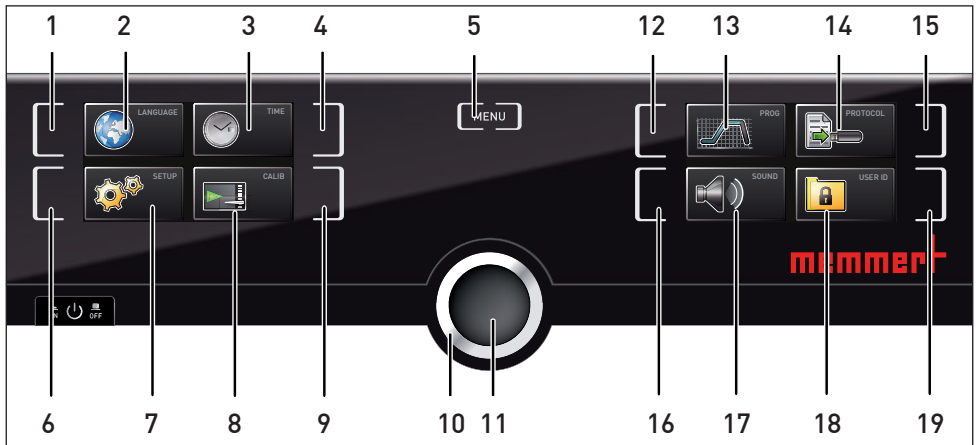



Fig. 24 ControlCOCKPIT in menu mode

- |  |  |
|--|--|
| <ul style="list-style-type: none"> <li>1 Language selection activation key</li> <li>2 Language selection display</li> <li>3 Date and time display</li> <li>4 Date and time setting activation key</li> <li>5 Exit menu mode and return to manual mode</li> <li>6 Setup activation key (basic appliance settings)</li> <li>7 Setup display (basic appliance settings)</li> <li>8 Adjustment display</li> <li>9 Adjustment activation key</li> <li>10 Turn control for adjustment</li> </ul> | <ul style="list-style-type: none"> <li>11 Confirmation key (accepts setting made with the turn control)</li> <li>12 Programme setup activation key</li> <li>13 Programme setup display</li> <li>14 Protocol display</li> <li>15 Protocol activation key</li> <li>16 Acoustic signal adjustment activation key</li> <li>17 Acoustic signal adjustment display</li> <li>18 User ID display</li> <li>19 User ID activation key</li> </ul> |
|--|--|

## 7.2 Basic operation in menu mode using the example of language selection

In general, all settings in menu mode are done just like in manual mode: Activate the respective display, use the turn control for setting and press the confirmation key to accept the change. A more detailed description is provided in the following, using the example of language selection.

1. Activate the desired parameter (in this example the language). To do so, press the corresponding activation key on the left or right of the respective display. The activated display is enlarged.

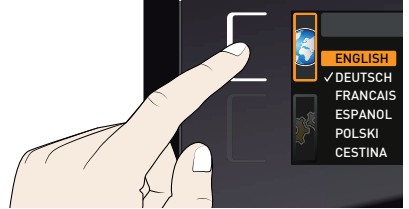



- 1 **i** If you want to exit or cancel your settings, again press the activation key which you have used to activate the display. The appliance returns to the menu overview. Only the settings that you have confirmed by pressing the confirmation key before cancelling the setting procedure are accepted.



2. With the turn control, select the desired new setting, e.g. Español (Spanish).



3. Save the setting by pressing the confirmation key.

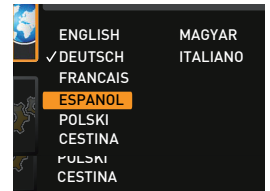
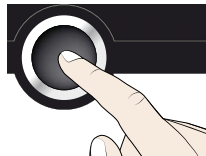


4. To return to the menu overview, press the activation key again.



You can now

- ▶ activate another menu function by pressing the corresponding activation key or
- ▶ return to manual mode by pressing the MENU key.

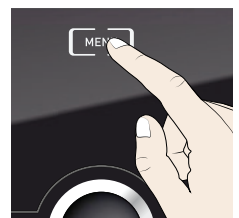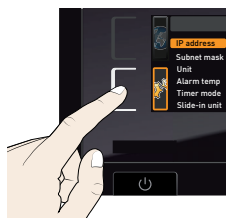

All other settings can be made accordingly. The settings possible are described in the following sections.

- 1 If no new values are entered or confirmed for approx. 30 seconds, the appliance automatically returns to the main menu and restores the former values.

## 7.3 Setup


### 7.3.1 Overview

In the SETUP menu, you can set the following parameters:

- ▶ the IP address and Subnet mask of the appliance's Ethernet interface (for connection to a network)
- ▶ the Unit of the temperature display (°C or °F, see page 45)
- ▶ Alarm temp: the temperature protection class according to DIN 12 880:2007-5 (TWW or TWB, see pages 45 and 33)
- ▶ how the digital backwards counter with target time setting works (Timer mode, see page 46)
- ▶ the type of the slide-in unit (Grid or Shelf, see page 44)
- ▶ the heat output distribution (Balance, see page 46)
- ▶ Remote control (see page 47)
- ▶ Gateway (see page 48)

- 1 If the SETUP menu contains more entries than can be displayed, this is indicated by the display "1/2". This means that there is a second „page“ of entries.

To display the hidden entries, use the turn control to scroll beyond the lowest entry. The page display changes to "2/2".

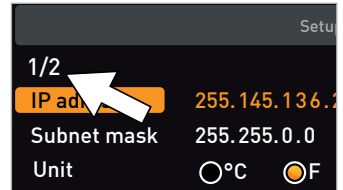

### 7.3.2 IP address and subnet mask

If you want to operate one or more appliances in a network, each appliance must have its own unique IP address for identification. By default, each appliance is delivered with the IP address 192.168.100.100.



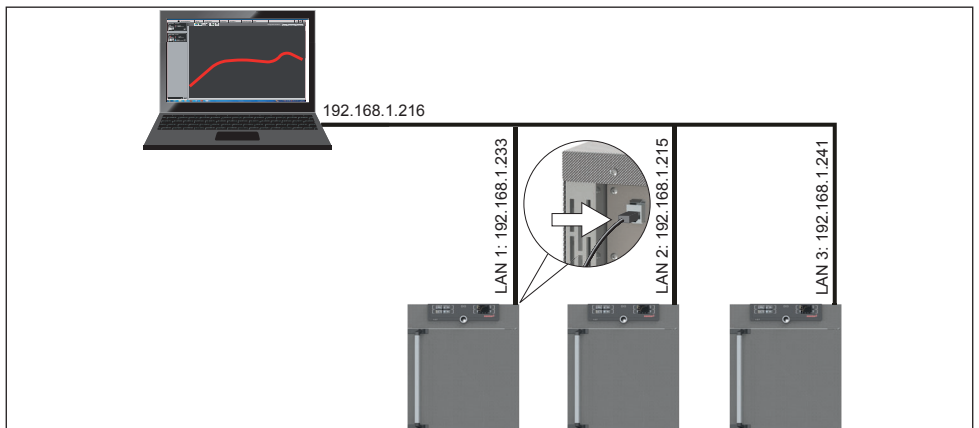


Fig. 25 Operation of several appliances in a network (schematic example)


1. Activate the SETUP display. The entry IP address is automatically highlighted.

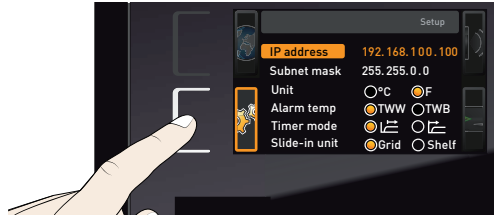


2. Accept the selection by pressing the confirmation key. The first three digits of the IP address are automatically selected.

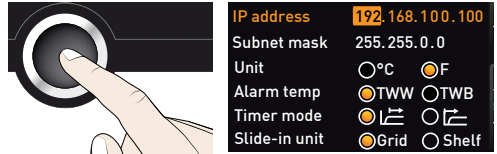


3. With the turn control, set the new number, e.g. 255.

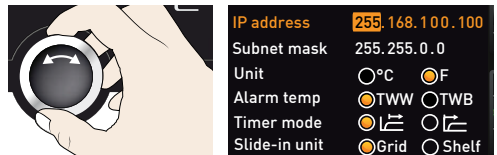

4. Accept the selection by pressing the confirmation key. The next three digits of the IP address are automatically selected. Setting these is done according to the description above.

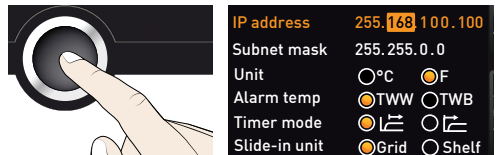

5. After setting the last three digits, accept the new IP address by pressing the confirmation key. The selection returns to the overview.

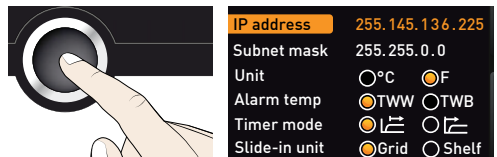

The subnet mask is set accordingly.

### 7.3.3 Unit

Here, you can choose whether the temperature is displayed in °C or °F.



### 7.3.4 Temperature monitoring (Alarm Temp)

Here, you can choose which temperature protection class in accordance with DIN 12 880:2007-5 should be used (TWW or TWB, description from page 33).



### 7.3.5 Timer mode

Here, you can choose whether the digital backwards counter with target time setting (see page 30, timer) should run setpoint-dependent or not – this determines whether the timer should not start until a tolerance band of  $\pm 3$  K around the set temperature is reached (Fig. 26, B) or whether it should start right after activation (A).

IP address	255.145.136.225	
Subnet mask	255.255.0.0	
Unit	<input checked="" type="radio"/> °C	<input type="radio"/> °F
Alarm temp	<input type="radio"/> TWW	<input checked="" type="radio"/> TWB
Timer mode	<input checked="" type="radio"/>	<input type="radio"/>
Slide-in unit	<input checked="" type="radio"/> Grid	<input type="radio"/> Shelf

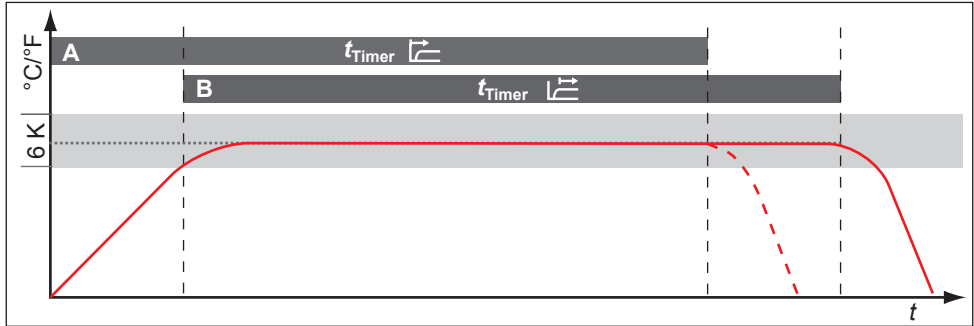


Fig. 26 Timer mode

- A Timer independent of setpoint: Timer starts right after activation
- B Timer setpoint-dependent: Timer does not start until tolerance band is reached

**i** If the temperature tolerance band is exceeded in setpoint-dependent operation, the timer will be interrupted and only be resumed when the setpoint temperature has been reached again.

### 7.3.6 Type of the slide-in unit (Grid or Shelf)

Here, you have to set the type of the slide-in unit (grid or shelf) used. The selection Shelf enables you to adjust the control function to the different air flow characteristics in the interior when using optional sliding shelves instead of the grids that are part of the standard delivery.

IP address	255.145.136.225	
Subnet mask	255.255.0.0	
Unit	<input checked="" type="radio"/> °C	<input type="radio"/> °F
Alarm temp	<input type="radio"/> TWW	<input checked="" type="radio"/> TWB
Timer mode	<input checked="" type="radio"/>	<input type="radio"/>
Slide-in unit	<input checked="" type="radio"/> Grid	<input type="radio"/> Shelf

### 7.3.7 Balance

For appliances of the size 55 and above, application-specific correction of the heat output distribution (balance) between the upper and lower heating groups is possible. The adjustment range is from  $-50\%$  to  $+50\%$ .

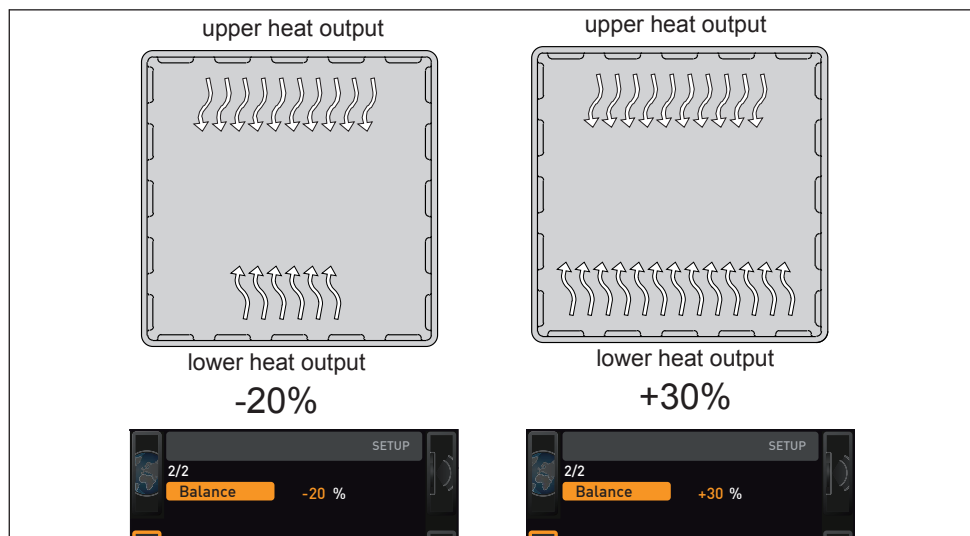




Fig. 27 Heat output distribution (example): The  $-20\%$  (left) setting causes the lower heating groups to emit  $20\%$  less heat than the upper ones. The  $+30\%$  (right) setting causes the lower heating groups to emit  $30\%$  more heat than the upper ones. The  $0\%$  setting restores the default heat output distribution.



### 7.3.8 Remote control

In the setup entry Remote control, you can set whether the appliance should be controlled via remote control and if so, in which mode. These settings are available:

- ▶ Off
- ▶ Read Only
- ▶ Write + Read
- ▶ Write + Alarm

When the appliance is in remote control mode, the  symbol appears in the temperature display. In the settings Write + Read and Write + Alarm, the appliance cannot be controlled at the ControlCOCKPIT until the remote control has been switched off (setting Off) or set to Read Only.


 In order to use the remote control function, programming skills and special libraries are required.

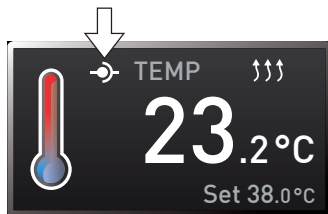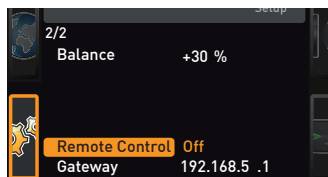

### 7.3.9 Gateway

The setup entry Gateway is used to connect two networks with different protocols.

The gateway is set the same way as the IP address (see page 44).

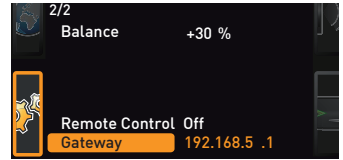


## 7.4 Date and Time


In the Time display, you can set date and time, time zone and daylight saving time. Changes can only be made in manual operating mode.

- 1 Always set the time zone (and summer time yes/no) before you set the date and time.
- 1 Avoid changing the set time after that since this can lead to gaps or overlapping when recording measured values. If you still need to change the time, you should not run a programme immediately before or after doing so.


1. Activate the time setting. To do so, press the activation key on the right side of the Time display. The display is enlarged and the first adjustment option (Date) automatically highlighted.

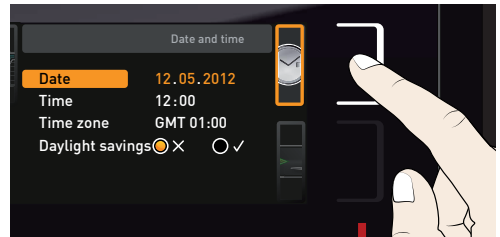


2. Turn the turn control until Time zone is highlighted.

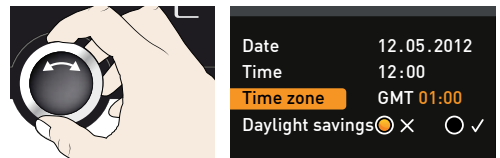


3. Accept the selection by pressing the confirmation key.

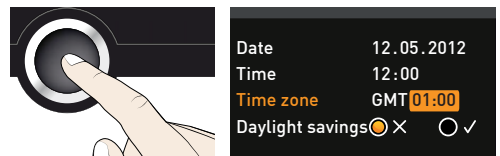

4. Set the time zone of the installation site with the turn control, e.g. 00:00 for Great Britain, 01:00 for France, Spain or Germany. Accept the selection by pressing the confirmation key.

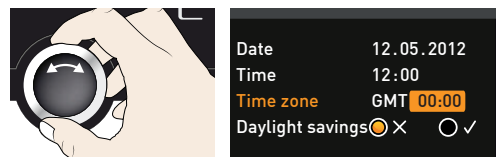


5. With the turn control, select the Daylight savings entry

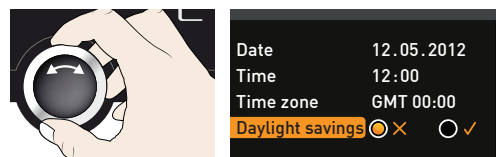




- Accept the selection by pressing the confirmation key. The adjustment options are highlighted.

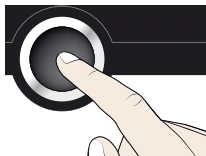


- Set daylight savings to off (X) or on (✓) with the turn control – in this case on (✓). Save the setting by pressing the confirmation key.

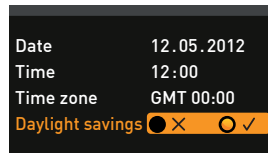

**i** Daylight saving time and standard time are not changed automatically. For this reason, please keep in mind to adjust them at the beginning of each period.

- Now, set date (day, month year) and time (hours, minutes) in the same way. Accept the selection by pressing the confirmation key.

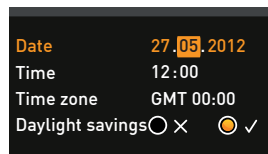

## 7.5 Calibration

The appliances are temperature calibrated and adjusted at the factory. In case readjustment should be necessary later on – for example due to influence of the chamber load – the appliance can be calibrated customer-specifically using three calibration temperatures of your choice:

- ▶ Cal1 Temperature calibration at low temperature
- ▶ Cal2 Temperature calibration at medium temperature
- ▶ Cal3 Temperature calibration at high temperature

To guarantee perfect control, we recommend to calibrate the appliance once a year.

**i** For temperature adjustment, you will need a calibrated reference measuring device.


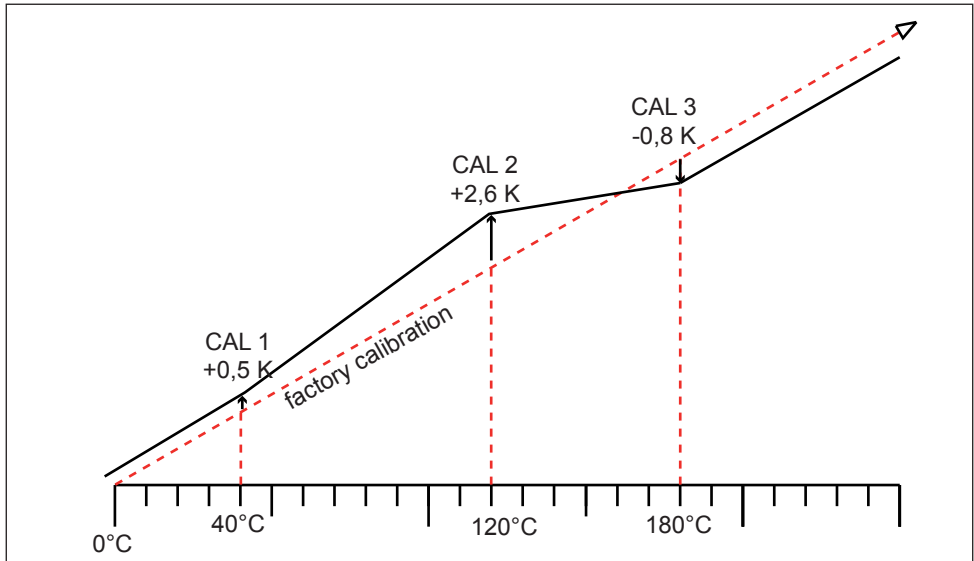



Fig. 28 Schematic example of temperature adjustment

Example: Temperature deviation at 120 °C should be corrected.

1. Activate the adjustment setting. To do so, press the activation key on the right side of the CALIB display. The display is enlarged and the first calibration temperature – in this case 40 °C – automatically highlighted.

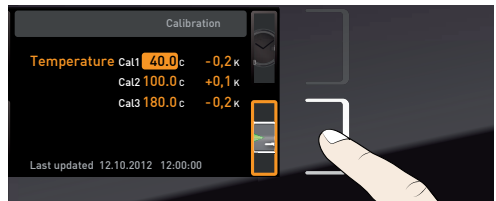

2. Press the confirmation key repeatedly, until the calibration temperature Cal2 is selected.

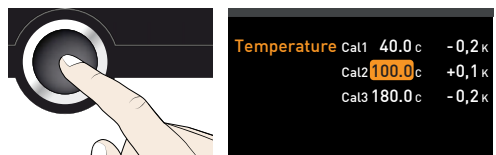


3. With the turn control, set the calibration temperature Cal2 to 120 °C.



Temperature	Cal1	40.0 c	-0,2 K
	Cal2	120.0 c	+0,1 K
	Cal3	180.0 c	-0,2 K

4. Save the setting by pressing the confirmation key. The corresponding calibration value is automatically highlighted.

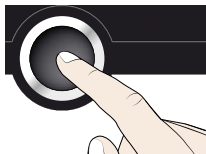

Temperature	Cal1	40.0 c	-0,2 K
	Cal2	120.0 c	+0,1 K
	Cal3	180.0 c	-0,2 K


5. Set the calibration value to 0.0 K and accept the setting by pressing the confirmation key.



Calibration			
Temperature	Cal1	40.0 c	-0,2 K
	Cal2	120.0 c	0,0 K
	Cal3	180.0 c	-0,2 K

6. Position the sensor of a calibrated reference instrument centrally in the appliance's working chamber.

7. Close the door and, in manual mode, adjust the set temperature to 120 °C.

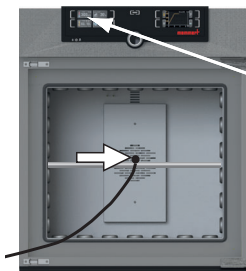

TEMP	21.4 °C
Set	120.0 °C

8. Wait until the appliance reaches the set temperature and displays 120 °C. The reference instrument for example displays 122.6 °C.

TEMP	120.0 °C
Set	120.0 °C

122.6 °C
----------

9. In the SETUP, adjust the calibration value Cal2 to +2.6 K (actual value measured minus setpoint temperature) and save the setting by pressing the confirmation key.



Temperature	Cal1	40.0 c	-0,2 K
	Cal2	120.0 c	+2,6 K
	Cal3	180.0 c	-0,2 K

10. After the calibration procedure, the temperature measured by the reference instrument should now also be 120 °C.

TEMP	120.0 °C
Set	120.0 °C

120.0 °C
----------


With Cal1, a calibration temperature below Cal2 can be programmed accordingly, and with Cal3, a temperature above. The minimum difference between the Cal values is 20 K for universal ovens UN../UF... and 10 K for incubators IN../IF...

**i** If all calibration values are set to 0.0 K, the factory calibration settings are restored.


## 7.6 Programme

In the PROG display, you can transfer programmes created using AtmoCONTROL software and saved on a USB storage medium to the appliance. Here, you can also select the programme to be used in manual mode (see page 31) and delete programmes.


**i** To load a programme from a USB storage medium : Connect the USB storage medium with the saved programme(s) to the interface on the right side of the control panel.

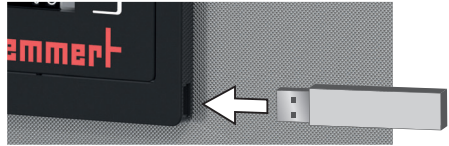


1. Activate the programme display. To do so, press the activation key on the left side of the PROG display. The display is enlarged and the entry Select automatically highlighted. The programmes available for activation are shown on the right. The programme currently available for use – in this example Test 012 – is highlighted in orange.

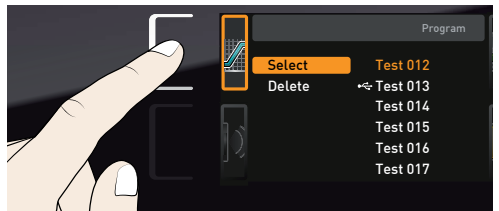


2. Call the Select function by pressing the confirmation key. All programmes available are displayed, including the ones saved on the USB storage medium (identified by the USB symbol). The programme currently available for use is highlighted in orange.

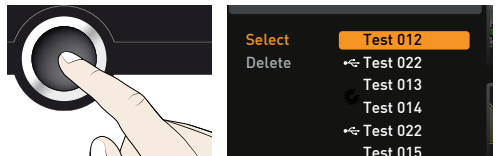


3. With the turn control, select the programme you want to make available for use.

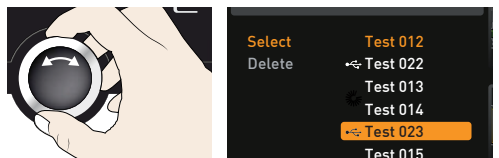

4. Accept the selection by pressing the confirmation key. The programme is now loaded, which is indicated by the transfer symbol.

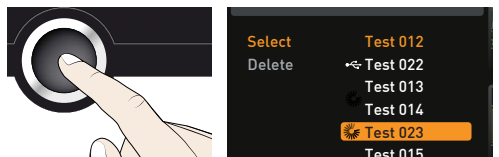

5. As soon as the programme is ready, the selection returns to Select. To start the programme: Return to manual mode by pressing the MENU key and start programme as described on page 31.

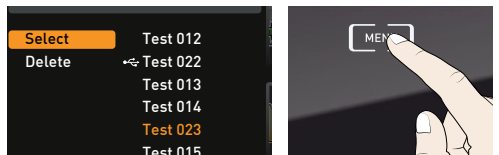

You can now remove the USB storage medium.

To delete a programme, select Delete with the turn control and select the programme to be deleted the same way you can select a programme for activation.

**i** A program to sterilise the work area in the controller is already pre-installed for Iplus appliances.


▶ Run time four hours at 160°C

## 7.7 Sound


In the SOUND display, you can define whether or not the appliance should emit acoustic signals and, if yes, define on which events it should do so:

- ▶ on the press of a key
- ▶ at the end of a programme
- ▶ on alarm
- ▶ if the door is open


1. Activate the acoustic signal adjustment. To do so, press the activation key on the left side of the SOUND display. The display is enlarged. The first category (in this case Key sound) is automatically highlighted. On the right, the current settings are shown on.

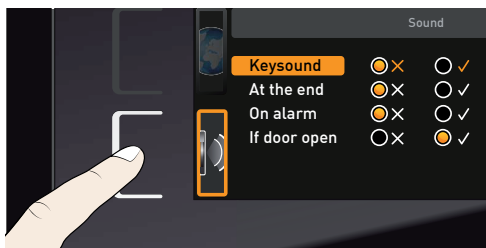


- i** If you want to edit another list entry: Turn the turn control until the respective entry – e.g. If door open (Special accessories) – is highlighted in colour.

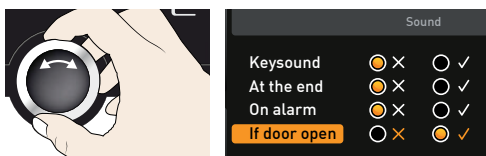

2. Save the selection by pressing the confirmation key. The adjustment options are automatically highlighted.

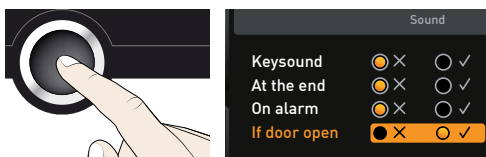


3. With the turn control, select the desired setting – in this example OFF (X).

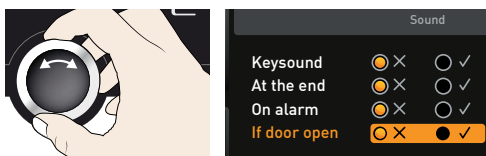

4. Save the setting by pressing the confirmation key.

- i** If an acoustic alarm sounds, it can be turned off by pressing the confirmation key.

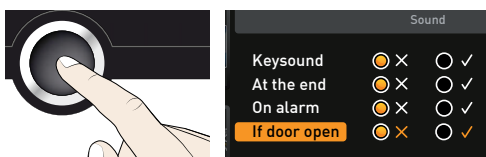


## 7.8 Protocol

The appliance continually logs all relevant measured values, settings and error messages at 1-minute intervals. The internal log memory is of the continuous memory type. The logging function cannot be switched off and is always active. The measured data are stored in the appliance, safe from manipulation. If the power supply is interrupted, the time of the power failure and voltage recovery are stored in the appliance.


You can export the protocol data for different periods to a USB storage medium via the USB interface or, via Ethernet, import them to the AtmoCONTROL software for graphical representation, print-out or storage.

**i** The log memory of the appliance is not modified or deleted by reading it out.


1. Connect the USB storage medium to the interface on the right side of the control panel.

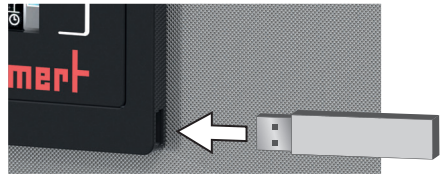


2. Activate the protocol. To do so, press the activation key on the right side of the PROTOCOL display. The display is enlarged and the period This Month automatically highlighted. To select another logging period, use the turn control.

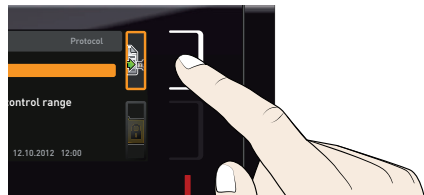

3. Save your selection by pressing the confirmation key. The transfer starts and a status symbol indicates the progress.

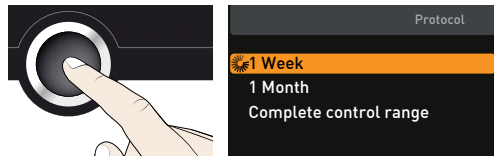

4. As soon as the transfer is complete, a check mark appears in front of the period selected. You can now remove the USB storage medium.

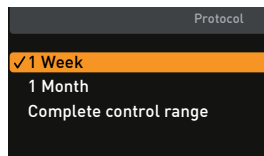

For a description of how to import and process protocol data in AtmoCONTROL or read them out via Ethernet, please observe the separate AtmoCONTROL manual.

## 7.9 User ID

### 7.9.1 Description

With the User-ID function, you can lock the settings of individual (e.g. temperature) or all parameters, so that they cannot be changed at the appliance by accident or unauthorised persons. You can also lock setting options in menu mode (e.g. adjustment or date and time settings) this way.

● If adjustment options are locked, this is indicated by **1** the lock symbol in the respective display (Fig. 29).

User ID data are entered in the AtmoCONTROL software and saved on the USB storage medium. The USB storage medium is thus acting as a key: Parameters can only be locked or unlocked if it is connected.

● A description of how to create a user ID in AtmoCONTROL is provided in the separate AtmoCONTROL manual.


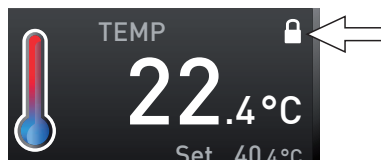



Fig. 29 Temperature adjustment at appliance locked (example)

### 7.9.2 User ID activation and deactivation


1. Connect the USB storage medium with the user ID data to the interface on the right side of the control panel.

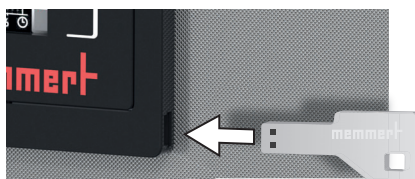

2. Activate the user ID. To do so, press the activation key on the right side of the USER-ID display. The display is enlarged and the entry Activate automatically highlighted.

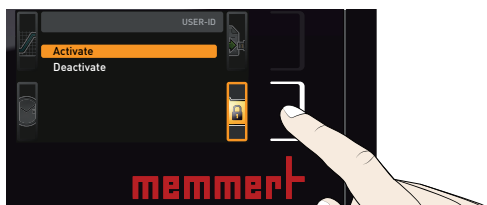

3. Confirm the activation by pressing the confirmation key. The new user ID data are transferred from the USB storage medium and activated. As soon as activation is complete, a check mark appears in front of the corresponding entry.

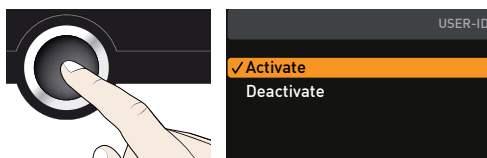

4. Remove the USB storage medium. Locked parameters are indicated by the lock symbol in the respective display (Fig. 29).

To unlock the appliance, connect the USB storage medium, activate the USER-ID entry and select Deactivate.

## 8. Maintenance and service

### 8.1 Cleaning



**Warning!**

**Danger due to electric shock. Before doing any maintenance work, pull out the mains plug.**



**Warning!**

**In case of appliances of a certain size, you can get accidentally locked in, which is life-threatening. Do not climb into the appliance!**



**Caution!**

**Danger of cuts due to sharp edges. Always wear gloves when working in the chamber interior.**

#### 8.1.1 Working chamber and metal surfaces

Regular cleaning of the easy-to-clean working chamber prevents build up of material remains that could impair the appearance and functionality of the stainless steel working chamber over time.

The metal surfaces of the appliance can be cleaned with normal stainless steel cleaning agents. Make sure that no rusty objects come into contact with the working chamber or with the stainless steel housing. Rust deposits can lead to an infection of the stainless steel. If rust spots should appear on the surface of the working chamber due to impurities, the affected area must be immediately cleaned and polished.

#### 8.1.2 Plastic parts

Do not clean the ControlCOCKPIT and other plastic parts of the appliance with caustic or solvent-based cleaning agents.

#### 8.1.3 Glass surfaces

Glass surfaces can be cleaned with a commercially available glass cleaner.

## 8.2 Regular maintenance

Once a year, grease the moving parts of the doors (hinges and lock) with thin silicone grease and check that the hinge screws are not loose.

To guarantee perfect control, we recommend to calibrate the appliance once a year (see page 50).

### 8.3 Repairs and service



**Warning!**

**After removing covers, live parts may be exposed. You may receive an electric shock if you touch these parts. Disconnect the mains plug before removing any covers. Any work inside the appliance may only be performed by qualified electricians.**



Repairs and service work are described in a separate service manual.



## 9. Storage and disposal

### 9.1 Storage

The appliance may only be stored under the following conditions:

- ▶ in a dry and enclosed, dust-free room
- ▶ frost-free
- ▶ disconnected from the power supply

### 9.2 Disposal

This product is subject to the Directive 2012/19/EC on Waste Electrical and Electronic Equipment (WEEE) of the European Parliament and of the Council. This appliance has been brought to market after August 13<sup>th</sup>, 2005 in countries which have already integrated this directive into their national laws. It may not be disposed of in normal household waste. For disposal, please contact your dealer or the manufacturer. Any appliances that are infected, infectious or contaminated with materials hazardous to health are excluded from return. Please also observe all other regulations applicable in this context.

Before disposing of the appliance, please render the door locking mechanism unusable, for example, to prevent playing children from being locked inside the appliance.

There is a lithium battery in the ControlCOCKPIT of the appliance. Remove it and dispose of it in accordance with the regulations in your country (Fig. 30).

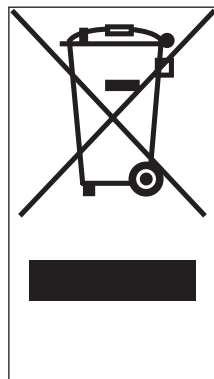


Fig. 30 Removing the lithium battery

#### Note for Germany:

The appliance may not be left at public or communal recycling or collection points.

## Index

- A**
- Accessories 16
  - Acoustic signals 50
  - Activation key 28
  - Adjustment 50
  - Air flap position 29
  - Air supply 11
  - Alarm 35, 38
  - Ambient conditions 15
  - Ambient temperature 15
  - Appliance error 39
  - ASF 33, 35
  - AtmoCONTROL 3, 12, 16, 28, 31, 52, 54, 55
  - Automatic temperature monitor 34
- B**
- Basic device settings 42
  - Basic settings 42
- C**
- CALIB 50
  - Calibration 50
  - Carrying 17
  - Cause of errors 39
  - Chamber load 26
  - Changes 8
  - Cleaning 56
  - Clock time 48
  - Communication interfaces 12
  - Compensation correction value 52
  - Connections 12
  - ControlCOCKPIT 9, 27
  - Convection 11
  - Customer service 2
- D**
- Declaration of conformity 15
  - Decommissioning 57
  - Delete programme 53
  - Delivery 17, 18, 23
  - Digital backwards counter 30
  - Dimensions 15
  - Disposal 57
  - Door 25
- E**
- Electrical connection 12
  - Electronic temperature monitoring 34
  - Emergency 8
  - Ending operation 37
  - End of programme 32
  - Error description 39
  - Error messages 38
  - Ethernet 12
- F**
- factory calibration settings 52
  - Fan speed 29
  - Fehlermeldung 40
  - Forklift truck 18
  - Function 10
- G**
- Gas bottles 23
  - Gateway 48
  - Graph 37
  - Grid 46
- H**
- Hazards 7
  - Heat output distribution 47
- I**
- Installation site 19
  - Intended use 10
  - Intended use as a medical device 10
  - Interfaces 12
  - Interior lighting 29
  - IP address 44
- L**
- Language selection 43
  - Lighting 29
  - Loading the appliance 26
  - Log memory 41, 54
- M**
- Maintenance 56
  - Malfunctions 8, 38
  - Manufacturer 2
  - Material 11
  - Mechanical temperature monitoring 35
  - Medical device 10
  - Menu 42, 48
  - Minimum clearances 19
  - Monitoring temperature 33
- N**
- Nameplate 13
  - Network 12, 44
  - Normal mode 28, 29
- O**
- Operating 25
  - Operating modes 28
  - Operating personnel 7, 25
  - Operating problems 39
  - Operation 25
- P**
- Packaging material 18
  - Power failure 41
  - Power supply 54
  - Press of a key 53
  - Product safety 7
  - Programme 52
  - Programme mode 28, 31
  - Pt100 temperature sensor 33
  - Putting into operation 23
- R**
- Rectifying errors 39
  - Regular maintenance 56
  - Remote control 47
- S**
- Safety regulations 6
  - Service 56
  - Servicing 56
  - Setting parameters 28, 43
  - Setting up 17, 19
  - Setting up options 20
  - Setup 44
  - Shelf 46
  - Slide-in unit 46
  - Speaker symbol 38
  - Storage after delivery 18
  - Switching off 37
  - Switching on 24

**T**

TB 35  
Technical data 14  
Temperature 29  
Temperature comparison 50  
Temperature deviation 50  
Temperature limiter 35  
Temperature monitor 33, 35  
Temperature monitoring 32,  
33, 45  
Temperature sensor 33  
Tilt protection 21  
Time 46  
Timer 30  
Timer mode 46  
Transport 17, 18  
Transport damage 18  
Türen einstellen 22  
Turn control 28  
TWB 34  
TWW 34  
TWW temperature monitor-  
ing 34

**U**

Unit 45  
Unpacking 18  
USB interface 12, 54  
User ID 55

**W**

Warning messages 12, 38  
Weight 14

# memmert

Universal ovens plus

Incubators plus

D33355 | Date 01/2020

englisch

Memmert GmbH + Co. KG  
Postfach 1720 | D-91107 Schwabach  
Tel. +49 9122 925-0 | Fax +49 9122 14585  
E-Mail: [sales@memmert.com](mailto:sales@memmert.com)  
[facebook.com/memmert.family](https://www.facebook.com/memmert.family)